


# The Instruction Set of the TANGO Controller

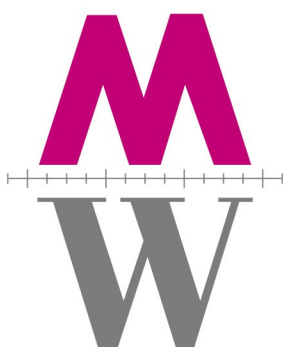

**MÄRZHÄUSER** WETZLAR

GMBH u. Co. KG

**Positioniersysteme**

In der Murch 15  
35579 Wetzlar  
Germany  
Tel.: +49/6441/9116-0  
[www.marzhauser.com](http://www.marzhauser.com)

# 1. Table of Contents

1. Table of Contents.....	2
2. Introduction.....	5
3. Brief Description of the Tango Instruction Set.....	6
4. Instruction Syntax Description.....	11
5. Error Numbers and their possible Root Cause.....	11
6. Controller Informations.....	12
6.1. version (Read detailed Version information).....	12
6.2. det (Read detailed Configuration).....	13
6.3. readsn (Read Serial Number).....	13
6.4. ver (Read default Version Number).....	14
6.5. iver (Read internal Version Number).....	14
7. Communication Interface Settings.....	15
7.1. baud (Baud Rate).....	15
7.2. cts (Enable/Disable RS232 Hardware Handshake).....	15
8. System Instructions.....	16
8.1. save (Save Parameters).....	16
8.2. restore (Restore Saved Parameters).....	16
8.3. reset (Force a Software Reset).....	16
8.4. pa (Enable or Disable the Power Amplifiers).....	17
8.5. ipreter (Select Instruction Set).....	17
9. Controller States and Error Messages.....	18
9.1. autostatus (Set Autostatus to required behavior).....	18
9.2. statusaxis (Query State of Axis).....	18
9.3. status (Query the Controller Error State).....	19
9.4. err (Query Error Number).....	19
9.5. help (Query Error Number with Description String).....	19
9.6. service (Print Service Information to Terminal).....	19
9.7. pci (Is PCI Bus).....	20
10. General Adjustments.....	21
10.1. dim (Unit for Positions).....	21
10.2. pitch (Spindle Pitch).....	21
10.3. gear (Gear Ratio).....	22
10.4. motorsteps (Motor Steps Per Revolution).....	22
10.5. accel (Maximum Acceleration).....	22
10.6. vel (Velocity).....	23
10.7. velfac (Velocity Factor).....	23
10.8. secvel (Secure Velocity).....	23
10.9. maxcur (Query Maximum Motor Current).....	24
10.10. cur (Motor Current).....	24
10.11. reduction (Motor Current Reduction Factor).....	24
10.12. curdelay (Delay for Current Reduction).....	25
10.13. axis (Enable, Disable, Switch Off Axis).....	25
10.14. axisdir (Axis Direction).....	26
10.15. stopaccel (Emergency Stop Deceleration).....	26
10.16. motortable (Motor Correction Table).....	26
11. Limit Switch Instructions (Hardware and Software).....	27
11.1. lim (Software Limits).....	27

11.2. limctr (Enable or Disable Limit Control).....	27
11.3. nosetlimit (Do not set limits by cal/rm).....	28
11.4. swtyp (Type of Limit Switch).....	28
11.5. swpol (Polarity of Limit Switch).....	29
11.6. swact (enable or disable limit switches).....	29
11.7. readsw (Read Status of Limit Switches).....	30
11.8. swin (Read Limit Switch Input Level).....	30
11.9. statuslimit (Limit Status).....	31
12. Calibration and Range Measure Instructions.....	32
12.1. cal (Command a Calibration).....	32
12.2. rm (Command a Range Measure).....	32
12.3. caliboffset (Calibration Offset).....	33
12.4. rmooffset (Range Measure Position Offset).....	33
12.5. caldir (Calibration Direction).....	33
12.6. calbspeed (Calibration Speed for Retraction).....	34
12.7. calrefspeed (Reference Signal Calibration Speed).....	34
12.8. calpos (Calibration Position).....	34
12.9. rekdir (Direction for Searching Reference Signal).....	35
13. Move Instructions.....	36
13.1. moa (Move Absolute).....	36
13.2. mor (Move Relative).....	36
13.3. m (Move Relative Shortcut).....	37
13.4. distance (Distance for m).....	37
13.5. moc (Move to Center).....	37
13.6. speed (Speed Move).....	38
13.7. a (Abort the Current Move).....	38
13.8. delay (Set the Delay Time for Consecutive Moves).....	38
13.9. pause (Set the Pause after Position Reached).....	39
13.10. pos (Read or Set Position).....	39
14. Joystick, Tackball and Handwheel Instructions.....	40
14.1. joy (Generally Enable/Disable/Set Joystick Mode).....	40
14.2. joydir (Joystick Direction or Assign Joystick).....	40
14.3. joywindow (Joystick Window).....	41
14.4. joyspeed (Joystick Speed Presets for BPZ Device).....	41
14.5. key (Read HDI Device Key State).....	41
14.6. hwfactor (Handwheel Transmission Factor).....	42
14.7. hwfilter (Handwheel Noise Filter).....	42
14.8. tbfactor (Handwheel Transmission Factor).....	42
14.9. tvrjoy (Pulse and Direction Joystick Functionality).....	43
14.10. tvrjoyf (Pulse and Direction Joystick Factor).....	43
15. Digital and Analogue I/O.....	44
15.1. digin (Digital Input).....	44
15.2. digout (Digital Output).....	44
15.3. anain (Analogue Input).....	45
15.4. anaout (Analogue Output).....	46
15.5. stoppol (Polarity of Stop Input Signal).....	46
15.6. shutter (Shutter Out Signal of AUX-IO).....	46
16. Encoder Instructions.....	47
16.1. encmask (Encoder Mask).....	47
16.2. enc (Encoder Active).....	47
16.3. encperiod (Encoder Signal Period).....	48



16.4. encdir (Encoder Counting Direction).....	48
16.5. encctl (Encoder has TTL Signal).....	48
16.6. encref (Use Encoder Reference Signal).....	49
16.7. encnas (Use Encoder NAS Error Signal).....	49
16.8. encnasstatus (Encoder NAS Error Signal State).....	50
16.9. encerr (Encoder Error State).....	50
16.10. encpos (Encoder Position).....	51
16.11. hwcount (Hardware Counter).....	51
16.12. clearhwcount (Clear Hardware Counter).....	51
17. MR Encoder Instructions.....	52
17.1. mra (MR Amplitude Correction Factor).....	52
17.2. mro (MR Amplitude Offset Correction Value).....	52
17.3. mrp (MR Signal Peak-To-Peak Measuring Result).....	53
17.4. mrt (MR Signal Level).....	53
18. Closed Loop Instructions.....	54
18.1. ctr (Control Enable).....	54
18.2. ctrc (Control Call).....	54
18.3. ctrt (Control Timeout).....	54
18.4. twi (Target Window).....	55
19. Trigger Signal Configuration.....	56
19.1. trig (Trigger).....	56
19.2. triga (Trigger Axis).....	56
19.3. trigm (Trigger Mode).....	57
19.4. trigger (Force Trigger Signal).....	58
19.5. trigs (Trigger Signal Length).....	58
19.6. trigd (Trigger Distance).....	58
19.7. trigf (Trigger Frequency).....	59
19.8. trigcount (Trigger Counter).....	59
20. Snapshot Signal Configuration.....	60
20.1. sns (Snapshot).....	60
20.2. snsl (Snapshot Level / Polarity).....	60
20.3. snsf (Snapshot Filter).....	61
20.4. snsm (Snapshot Mode).....	61
20.5. snsc (Snapshot Counter).....	61
20.6. snsp (Snapshot Position).....	62
20.7. snsarray (Snapshot Array).....	62
20.8. prehome (Snapshot PreHome Position).....	63
20.9. home (Snapshot Home Position).....	63
21. Scan Mode.....	64
21.1. scanmode (Switch to Scan Mode).....	64
21.2. scanvel (Scanmode Vector Velocity).....	64
22. Document Revision History.....	65

## 2. Introduction

All instructions and parameters which are sent to the controller, as well as all feedbacks of the controller, are transferred as a sequence of ASCII characters. The connection may be controlled manually at any time with a terminal program (e.g. HyperTerminal). The use of ASCII string communication simplifies error tracing, if the instructions are given over a customized program. Commands from the PC to the controller are marked with either an exclamation mark '!', if they transmit parameters to the controller or with a preceding question mark '?', if the PC requests data from the controller. The controller does not distinguish between upper and lower case ASCII characters. All floating point decimal numbers contain a point and no comma between pre and postpositions.

Examples:

```
!cal      command the controller to do a calibration
?status  ask for the status of the controller
```

### Hints:

In some cases, like single character instructions, it is not necessary to use the leading indicator '!' or '?'. Other instructions, e.g. for moving the axes, require the delivery of parameters. These are transferred following the instruction. Blanks must be inserted between the command text, its parameters and to separate each parameter. Example: "moa 45 13 20[CR]" means proceed x, y and z to the positions 45, 13 and 20. Each instruction must be terminated with a carriage return [CR]. This indication is represented as follows in the ASCII character set:

Symbol	decimal value	hexadecimal value	binary value
CR	13	0xD	00001101

Move commands are executed as a vector move, so they complete at the same time. To move axes independently, with their own velocities, they may be started by separate single axis commands. Please refer to the "move" command description.

Many parameters can be stored permanently in the Tango Controller, so they are available after each consecutive power on. When stored once, this reduces initialization overhead of the application software. Refer to the "save" command for further information. Parameters that are saved can be identified by a 'Y' in the Save column of the brief instruction description.

Please do not send more than 255 characters at once to the Tango Controller, as the input buffer will overflow. To avoid this it is recommended to request the "?err" state inbetween and wait for a value to be returned. Another solution is to activate the "!cts" handshake (available in Desktop RS232 or USB versions). This will automatically halt the PC transmission for as long as the input buffer is full. The PC COM port then must be opened with hardware handshake on, too. Please refer to the "!cts" command description.

### Important: Security speed limitation!

The Tango controller has a built in security function, which reduces the maximum travel velocity to 10mm/s for as long as no initial cal/rm move is executed. This is to preserve the microscope stage from damage that could be caused by moving fast into its end positions. After calibrating the axis into its endswitches (cal and/or rm if switches are mounted and enabled) the travel velocity is no longer limited.

If it is not wanted or impossible to do a calibration/range measure move on each power on, the speed limit may be increased to up to 100mm/s at own risk. Please refer to the "secvel" command for further information.

### 3. Brief Description of the Tango Instruction Set

Controller Informations					
Instruction	Example	Save	Brief description	Page	
(?)	version	version	fix	Read detailed firmware and controller version	12
(?)	det	det	fix	Read detailed configuration information	13
?	readsn	?readsn	fix	Read the controller serial number	13
(?)	ver	ver	fix	Read default version number	14
(?)	iver	iver	fix	Read further version number information	14

Communication Interface Settings					
Instruction	Example	Save	Brief description	Page	
?!	baud	!baud 9600	Y	Set RS232 baud rate to 9600 Bd (default=57600)	15
?!	cts	!cts 1	Y	Switch on CTS hardware handshake	15

System Instructions					
Instruction	Example	Save	Brief description	Page	
(!)	save	save	-	Save present parameters to controller nonvolatile memory	16
(!)	restore	restore	-	Reset present parameters to saved values	16
(!)	reset	reset	-	Reset controller (forces restart, similar to cycle power)	16
!	pa	!pa 1	-	Enable power amplifiers (disable = 0), see 'axis' cmd. too	17
?!	ipreter	!ipreter 1	Y	Select standard LSTEP command set	17

Controller States and Error Messages					
Instruction	Example	Save	Brief description	Page	
?!	autostatus	!autostatus 0	-	Select autostatus response type 0 (=disabled), range: [04]	18
(?)	statusaxis	statusaxis	-	Read axis state [@,M,J,C,S,A,D,]	18
(?)	status	status	-	Read controller error state	19
(?)	err	err	-	Read error number	19
(?)	help	help	-	Read error number with additional text description	19
(?)	service	service	-	Returns a detailed parameter and state list, for debugging	19
(?)	pci	pci	-	Returns 1 if controller is plugged in a PCI slot (desktop=0)	20

General Adjustments					
Instruction	Example	Save	Brief description	Page	
?!	dim	!dim 1 1 1	Y	Set position units of X Y Z to $\mu\text{m}$	21
?!	pitch	!pitch 1 1 1	Y	Set spindle pitch of X Y Z to 1 [mm/revolution]	21
?!	gear	!gear 1 1 1	Y	Set gear factor of X Y Z to 1	22
?!	motorsteps	!motorsteps x 200	Y	Set X axis motor has 200s steps per revolution	22
?!	accel	!accel 0.1 0.1 0.1	Y	Set acceleration of X Y Z to $0.1\text{m/s}^2$	22
?!	vel	!vel 10 10 10	Y	Adjust speed of X Y Z to 10 [revolutions/s]	23
?!	velfac	!velfac 1 1 1	Y	Set velocity reduction factor for X Y Z to 1 (= no reduction), range is [0.01-1]	23
?!	secvel	!secvel x 20	Y	Set secure speed limit X to 20mm/s (unit is always mm/s)	23
?	maxcur	?maxcur	fix	Show the maximum possible motor currents of all axes	24
?!	cur	!cur 0.5 0.6 1	Y	Set motor current in Ampere: X=0.5 Y=0.6 and Z=1 A	24
?!	reduction	!reduction 0.5 0.5 0.5	Y	Select 50% motor current reduction for X Y Z	24
?!	curdelay	!curdelay 1000	Y	Delay X axis motor current reduction by 1000 [ms]	25
?!	axis	!axis 1 0 -1	Y	Enable X, disable Y and switch off Z axis	25
?!	axisdir	!axisdir 0 1 0	Y	Reverse rotating direction of Y motor (caution!)	26
?!	stopaccel	!stopaccel 2	Y	Set deceleration during stop condition to $2\text{m/s}^2$	26
?!	motortable	!motortable x 2	Y	Select custom motor correction table type 2 for X axis	26

Limit Switch Instructions (Hardware and Software)					
Instruction	Example	Save	Brief description	Page	
?!	lim	!lim 0 10 0 10 0 10	-	Set lower position limit to 0 and upper limit to 10 (assume unit is [mm] if dim was set to 2) for X Y Z	27
?!	limctr	!limctr x 1	-	Enable hardware limit switches for X axis, default = 1	27
?!	nosetlimit	!nosetlimit 1 1 1 1	Y	Disable setting/overwriting of software limits during cal and rm for all axes (here: X Y Z A), default = 0	28
?!	swtyp	!swtyp 1 0 1 !swtyp y 0 0 0	Y	Set limit switch type for all axes to NPN (pull-up) Set limit switch type for Y to PNP (pull-down)	28
?!	swpol	!swpol 1 0 1 !swpol z 1 0 1	Y	Set polarity of limit switches for all axes to active high (=1) Set polarity of limit switches for Z to active high	29
?!	swact	!swact 1 0 1 !swact y 1 0 0	Y	Enable cal and rm limit switches for all axes Enable cal limit switch for Y, disable ref and rm	29
?	readsw	?readsw	-	Read states of all limit switches (1=active and actuated)	30
(?)	swin	swin	-	Read TTL signal level of all limit switch inputs (1=high)	30
(?)	statuslimit	statuslimit	-	Read current limit status „A“ = calibration done „D“ = rm done „L“ = limit switch modified by software „“ = not yet modified	31

Calibration and Range Measure Instructions					
Instruction	Example	Save	Brief description	Page	
(!)	cal	cal	-	Perform a calibration move for all enabled axes, see 'axis'	32
(!)	rm	rm x	-	Perform a range measure move in X	32
?!	caliboffset	!caliboffset 1 1 1	Y	Set the cal zero-point 1mm aside lower limit switch (dim 2)	33
?!	rmoffset	!rmoffset 1 1 1	Y	Set rm end-position 1mm aside upper limit switch (dim 2)	33
?!	caldir	!caldir z 1	Y	Calibrate the Z-axis in positive direction	33
?!	calbspeed	!calbspeed 20	Y	Set the speed for move out of 'cal' and 'rm' limit switches for all axes to 0.2 [revolutions/s], range is [1...100]	34
?!	calrefspeed	!calrefspeed 10	Y	Set the speed for calibrating to the encoder reference for all axes to 0.1 [revolutions/s], range is [1...100]	34
?!	calpos	calpos	-	Read back the encoder position where the calibration switch was released	34
?!	refdir	?refdir y	-	Read the direction for encoder reference search in Y axis	35

Move Instructions					
Instruction	Example	Save	Brief description	Page	
(!)	moa	moa 10 10 10 moa y 20	-	Move X Y Z absolute to positions 10 10 10 Move Y axis to position 20 (unit depends on dim setting)	36
(!)	mor	mor 4 4 4 mor y -10.5	-	Move X Y Z relative by 4 (unit depends on dim setting) Move Y axis relative 10.5 backwards	36
(!)	m	m	-	Move relative again (use same parameters as defined by last '!mor' or '!distance' instruction)	37
?!	distance	!distance 1 1 1	-	Set distance for X Y Z 'm'-move (start with 'm' or '!m')	37
(!)	moc	moc x	-	Move X to center position between lower and upper limit switch, or between lower and upper software limits	37
?!	speed	!speed 5 5 5 !speed y 0	-	Digital joystick: move X Y Z axis with 5 [revolutions/s] Stop the Y axis speed move	38
(!)	a	a	-	Abort move (Stop)	38
?!	delay	!delay 1000	Y	Delay all consecutive moves by 1000 ms	38
?!	pause	!pause 10	Y	Delay "position reached" autostatus response by 10 ms	39



? !	pos	!pos 0 0 0 !pos z 1.2	-	Set current X Y Z position to 0 Set current Z position to 1.2	39
-----	-----	--------------------------	---	--	----



### Joystick, Tackball and Handwheel Instructions

Instruction	Example	Save	Brief description	Page	
?!	joy	!joy 0 !joy 2	Y	Switch joystick on(=2) or off(=0)	40
?!	joydir	!joydir 1 1 1	Y	Set motor direction for joystick operation (Z reversed)	40
?!	joywindow	!joywindow 14	Y	Set idle window of the joystick center position, where a joystick deflection has no effect [0..100]	41
? (!)	joyspeed	joyspeed 2 25	Y	Set joystick speed for speed button 2 "medium" to 25 rev/s	41
(?)	key	key	-	Read state of all joystick buttons (0=released, 1=pressed)	41
?!	hwfactor	!hwfactor x 100	Y	One handwheel revolution in X is 100mm stage travel	42
?!	hwfilter	!hwfilter 0	Y	Switch off handwheel noise reduction	42
?!	tbfactor	!tbfactor 1 1	Y	Set trackball transmission factor in X and Y to default	42
?!	tvrjoy	!tvrjoy z	Y	Assign AUX-IO pulse&direction joystick to Z axis	43
?!	tvrjoyf	!tvrjoyf 1	Y	Set tvrjoy transmission factor to 1	43

### Digital and Analogue I/O

Instruction	Example	Save	Brief description	Page	
(?)	digin	digin digin 8	-	Read all digital inputs Read digital input 8	44
?!	digout	!digout 5 1 ?digout	-	Set digital output 5 to logic level 1 Read back all digital output levels	44
(?)	anain	anain c 2	-	Read input of analogue channel 2	45
?!	anaout	!anaout c 1 17.5	-	Set analogue voltage of channel 1 to 17.5 percent (1.75V)	45
?!	stoppol	!stoppol 1	Y	AUX-IO stop input is active high (=1) or active low (=0)	46
?!	shutter	!shutter 1	-	Set AUX-IO shutter out signal to TTL high	46

### Encoder Instructions

Instruction	Example	Save	Brief description	Page	
?!	encmask	!encmask 1 1 0	Y	Enable activation of X and Y encoders, disable Z	47
?!	enc	!enc 1 0	-	Manually activate X encoder (caution!), set Y to inactive	47
?!	encperiod	!encperiod 0.1	Y	Set signal period of X encoder to 100 µm	48
?!	encttl	!encttl x 1	Y	X encoder is TTL type (has no analogue sin/cos signal)	48
?!	encdir	!encdir y 1	(Y)	Reverse counting direction for Y encoder	48
?!	encref	!encref 0	Y	No decoding of X encoder reference signal	49
?!	encnas	!encnas 1 0 0	Y	Enable NAS error signal input encoding for X encoder only	49
(?)	encnasstatus	encnasstatus x	-	Read X encoder NAS signal state (1=NAS error)	50
?!	encerr	!encerr 0	-	Clear encoder error state for X axis (? response is 0 or e)	50
?!	encpos	!encpos 1	-	?pos instruction returns for X the encoder position, if enc=1	51
(?)	hwcount	hwcount	-	Read all encoder positions (TTL counter, no interpolation)	51
(!)	clearhwcount	clearhwcount x	-	Set X axis hwcount to zero	51

### MR Encoder Instructions

Instruction	Example	Save	Brief description	Page	
?!	mra	?mra x	-	Read amplitude correction factor (sin/cos ratio) of X	52
?!	mro	?mro	-	Read offset correction value for all encoders	52
?!	mrp	!mrp x 0 0 0 0	-	Reset MR-signal peak-to-peak measurement result of X	53
?	mrt	?mrt z 2	-	List two measurement results of the Z input signals	53

Closed Loop Instructions					
Instruction	Example	Save	Brief description	Page	
?!	ctr	!ctr 1 1 1	Y	Set closed loop circuit to "active until reached" mode	54
?!	ctrc	!ctrc 1	Y	Closed loop control is called every 1 millisecond	54
?!	ctrt	!ctrt 5	Y	Closed loop control Timeout after 5 milliseconds	54
?!	twi	!twi 0.01 0.01 0.01	Y	Set target window for X Y Z to 10µm (assume dim=2)	55

Trigger Signal Configuration <sup>1</sup>					
Instruction	Example	Save	Brief description	Page	
?!	trig	!trig 1	-	Enable trigger functionality (should be the last command)	56
?!	triga	!triga x	-	Trigger function is related to X axis	56
?!	trigm	!trigm 0	-	Select trigger mode 0	57
(!)	trigger	trigger	-	Manually set trigger output (available in trigm 102, 103)	58
?!	trigs	!trigs 40	-	Set trigger output signal length to 40 microseconds	58
?!	trigd	!trigd 10	-	Set trigger distance to 10 (mm if dim=2)	58
?!	trigf	!trigf 1000	-	Generate periodic trigger pulses with 1kHz	59
?!	trigcount	?trigcount	-	Read number of generated trigger events	59

Snapshot Signal Configuration <sup>1</sup>					
Instruction	Example	Save	Brief description	Page	
?!	sns	!sns 1	-	Enable snapshot functionality (should be last command)	60
?!	snsi	!snsi 0	-	Set snapshot input signal to active low	60
?!	snsf	!snsf 10	-	Set snapshot signal glitch filter to 10 milliseconds	61
?!	snsm	!snsm 0	-	Set snapshot mode to 0(=capture, 1=move)	61
?!	snsr	?snsr	-	Read number of snapshot events (=array fill size)	61
?!	snsx	?snsx	-	Read last captured X position	62
?!	snsa	?snsa 1	-	Read first position entry of snapshot array (all axes)	62
?!	prehome	!prehome 10 20 1	-	Set prehome positions X Y Z to 10 20 1 (unit depends on dim setting)	63
?!	home	!home 5 5 0	-	Set home positions X Y Z to 5 5 0 (unit depends on dim setting)	63

Scan Mode					
Instruction	Example	Save	Brief description	Page	
?!	scanmode	!scanmode 1	-	Set controller behavior to scanmode	64
?!	scanvel	!scanvel 20 20	-	Set scanmode vector velocity to 20mm/s for X and Y	64

<sup>1</sup> Function has to be enabled by factory, it is not available per default.

## 4. Instruction Syntax Description

Most instructions work in both directions (reading and writing). (?)! means the instruction accepts write and read. The controller identifies a read command by the preceded '?', or '!' for writing parameters.

Some examples for legal instruction syntax:

```
(?)!Command parameter1 parameter2 parameter3 parameter4  
(?)!Command parameter1 parameter2  
(?)!Command axis parameter  
(?)!Command
```

## 5. Error Numbers and their possible Root Cause

```
0 no error  
1 no valid axis name  
2 no executable instruction  
3 too many characters in command line  
4 invalid instruction  
5 number is not inside allowed range  
6 wrong number of parameters  
7 either ! or ? is missing  
8 no TVR possible, while axis active  
9 no ON or OFF of axis possible, while TVR active  
10 function not configured  
11 no move instruction possible, while joystick enabled  
12 limit switch active  
13 function not executable, because encoder detected  
21 multiple axis moves are forbidden (e.g. during initialization)  
22 automatic or manual move is not allowed (e.g. door open or initialization)  
27 emergency STOP is active  
29 servo amplifier are disabled (switched OFF)  
30 safety circuit out of order
```

## 6. Controller Informations

You may read the firmware version by sending the instruction 'version' to the controller. The instruction 'det' gives you further details of which options are enabled. Each controller has its own unique serial number readable with the instruction 'readsn'.

### 6.1. version (Read detailed Version information)

Syntax: ?version or version

Parameter: none

Description: This instruction delivers detailed information about the firmware version.

Example: ?version

TANGO-DT-S, Version 1.24, Jun 20 2007 , 11:19:32

Response syntax: Character string including controller type, firmware version and build date separated by a comma:

TANGO Fixed string identifying the Tango controller

-DT Desktop version

-PCI PCI card version

-S Tango short card version (PCI-S, DT-S)

Version 1.24 Firmware version number

Jun 20 2007 Firmware build date

11:19:32 Firmware build time

## 6.2. det (Read detailed Configuration)

Syntax:           ?det or det  
 Parameter:       none  
 Description:      This instruction delivers detailed information about the current controller configuration.  
 Response:        The controller returns a decimal integer number. Its hexadecimal value represents the configuration, like following table shows:

```

0x0    1    1Vssencoder is configured
0x0    2    MREncoder is configured
0x0    4    TTLencoder is configured
0x0    3    this is the number of configured axes (e.g. 3)
0x0    1    Display is configured
0x0    2    Speedpoti is configured
0x0    4    Hand wheel is configured
0x0    8    Snapshot is configured
0x0    1    TVRin is configured
0x0    2    Trigger out is configured
0x0    8    TVRout is configured
0x1    16   digital I/O are configured
0x2    32   digital I/O are configured
0x4    1    Trackball is configured
  
```

The current configuration results as a logical 'or' of these bits.

Example:           Assume ?det delivers the response 81697 which is 13F21 hexadecimal. This number means in detail, that the controller is configured for:

```

1 => 16 digital I/O
3 => TVRin and Trigger out
F => Display, Speedpoti, Hand wheel and Snapshot
2 => 2 axis
1 => 1Vss encoder
  
```

## 6.3. readsn (Read Serial Number)

Syntax:           ?readsn or readsn  
 Parameter:        none  
 Description:      The instruction ?readsn delivers the current serial number.  
 Example:          ?readsn  
 Response:        The controller transmits its unique serial number as ASCII characters like YYWWNNXXX.

```

YY    year of manufacturing
DD    week of manufacturing
NN    available axes (in hardware)
XXX   Index number
  
```

## 6.4. ver (Read default Version Number)

Syntax: ?ver or ver  
Parameter: none

Description: This instruction reads back the default firmware version info. The first number is the number of configured axes. The second number is the maximum possible motor current in ampere. For Tango firmware version information please use "version".

Example: ?ver  
Response syntax: Vers:LSnm.xx.xxx

"Vers:LS"	Fixed character string
n	Number of configured axes: 1, 2, 3, or 4
m	Maximum Current: 1=1.25A, 2=2.5A, 3=3.75A
x	Fixed number

## 6.5. iver (Read internal Version Number)

Syntax: ?iver or iver  
Parameter: none

Description: This instruction reads the internal version information string. Mostly unused. Please use the "version" command to read the Tango firmware version.

Response syntax: 'T''day of week''.''week''.''year''-''number'  
Example of ?iver response: T04.35.020004

## 7. Communication Interface Settings

### 7.1. baud (Baud Rate)

Syntax:           !baud or ?baud  
Parameter:       9600, 19200, 38400, 57600 or 115200

Description:       This instruction sets or reads the serial communication transfer rate (baudrate). After sending this command first make sure the controlling device (e.g. a PC) has the same setting again. Then a save command may be sent to permanently store the new baudrate.  
For PCI bus communication this instruction has no effect.

Response:         Current baud rate.

Examples:  
!baud 57600        The baud rate is set to 57600 [Bd].  
?baud             query controller for current baud rate

### 7.2. cts (Enable/Disable RS232 Hardware Handshake)

Syntax:           ?cts or !cts  
Parameter:        0 or 1

Description:       Writing a 1 enables additional hardware handshake of the RS232 or USB interface. A 0 disables this function.  
For PCI bus communication this instruction has no effect.  
Please note that the PC COM port has to be opened in hardware handshake mode, too.

Response:         Current state of CTS (0=disabled or 1=enabled)

Examples:  
?cts              query controller for current state of CTS  
!cts 0            disable CTS handshake  
!cts 1            enable CTS handshake

## 8. System Instructions

The controller provides two different instruction sets.

> The default instruction set as described in this manual.

> The second optional instruction set is a subset of the Venus command set. The following instruction let you select your required instruction set (if the option is installed).

### 8.1. save (Save Parameters)

Syntax:               !save or save  
Parameter:           none

Description:        The instruction !save stores your favorite parameter settings (like spindle pitch) in a permanent and safe data area. These parameters will be taken by the controller after each consecutive reset or power on as default values. Executing a save comand always returns the "OK..." string when writing is completed.

Response:            The save instruction returns the response string "OK..."

Example:

save => The currently used controller parameters are saved as default.

### 8.2. restore (Restore Saved Parameters)

Syntax:               !restore or restore  
Parameter:           none

Description:        The controller reloads the saved parameters from its nonvolatile memory. The current controller parameters get overwritten by the saved defaults. Refer to the "save" instruction. Similar to a software "reset", but without affecting the hardware.

Response:            none  
Example:              restore

### 8.3. reset (Force a Software Reset)

There are two possibilities to reset the controller:

- . The power on reset
- . The Software Reset

Syntax:               !reset or reset  
Parameter:           none

Description:        The controller is forced to perform a software reset. It is a restart similar to power on. Rebooting from reset will take more than 1 second, where the controller is not responding. There is no reply to a software reset. So for knowing if the controller is rebooted and ready, it may be necessary to poll data until it responds again.

Response:            none  
Example:              reset



## 8.4. pa (Enable or Disable the Power Amplifiers)

Syntax:           !poweramplifier or !pa  
Parameter:       0 or 1

Description:      This instruction switches all motor amplifiers on(=1) or off(=0). If switched off, no motor current is flowing. To switch off axes individually, please use the 'axis' command.

Response:        none

Example:          !pa 1            Switch on all amplifiers.

## 8.5. ipreter (Select Instruction Set)

Syntax:           !ipreter or ?ipreter  
Parameter:        0, 1 or 2

Description:      0 => Prohibited. Register command set is no longer provided.  
                  1 => Default instruction set (LSTEP), as described in this manual.  
                  2 => Optional instruction set (VENUS-1).

To return from the VENUS instruction set (2), please enter the string "1 setipreter" and press enter (or send an ASCII [CR]).

Response:        0, 1 or 2 (here certainly 1 only)

Example:  
!ipreter 1 => The controller uses the LSTEP interpreter.  
?ipreter   => Responds the currently selected interpreter.

## 9. Controller States and Error Messages

### 9.1. autostatus (Set Autostatus to required behavior)

Syntax: !autostatus or ?autostatus

Parameter: 0, 1, 2, 3 or 4

Description:

- 0 => Sending of any automatic state messages is switched OFF, except 'save'.
- 1 => After each automatic move the message 'position reached' (this is the character '@' for each configured axis) is automatically transmitted by the controller. It is the default configuration after power on.
- 2 => The controller transmits the message 'position reached' plus other status messages.
- 3 => Instead of 'position reached' string a simple <CR> is transmitted (=fast).
- 4 => Echoes all input instructions including parameters.

Example: Assume there are three axes configured and autostatus is set to 1.  
After completion of a move (moa, mor, m, a) the controller will return a "@@@-." which means position reached.

```
!autostatus 0      Switch off autostatus (now the position reached info has to be
                  polled by using the "statusaxis" instruction).
?autostatus       Reads back the selected autostatus value.
```

### 9.2. statusaxis (Query State of Axis)

Syntax: ?statusaxis or statusaxis

Parameter: none

Description:

Statusaxis responds the state of each axis.  
Similar to the 'autostatus 1' response to move commands,  
but with an additional '-' after the dot.  
Can be used for polling axes in 'autostatus 0' mode.

```
@ => Axis is not moving and ready
M => Axis is moving
J => Axis is controlled manually (by joystick)
C => Axis is in closed loop
S => Limit switches are triggered and prevent further automatic move
A => ok response after cal instruction
D => ok response after rm instruction
E => not o.k. response after cal or rm, if an error occurred during cal
    instruction (e.g. a limitswitch is not working proper)
U => manual adjustment (e.g. 1st setup)
T => Timeout
=> Axis is not enabled
```

Example: Assume ?statusaxis delivers the response @@@-.-  
This means three axis enabled and read to move.

### 9.3. status (Query the Controller Error State)

Syntax:            ?status or status  
Parameter:        none

Description:       The ?status instruction responds with the current state of the controller. Which is either 'OK...' or an 'ERR' with error number. Also see 'err' instruction.

Response:         OK... or ERR with error number

Example:           ?status => ERR 4

### 9.4. err (Query Error Number)

Syntax:            ?err or err  
Parameter:        none

Description:       The instruction err returns the current error number or 0, if no error occurred.

Response:         Error number as decimal value  
                  (refer to Chapter 5 "Error Numbers")

Example:           err => 0

### 9.5. help (Query Error Number with Description String)

Syntax:            ?help or help  
Parameter:        none

Description:       The instruction help returns a text string. It contains the error state (current error number) like the 'err' instruction, but with an additional error description string. May be used for debugging with a terminal program.

Response:         Error number as decimal value, error description

Example:           help => ERROR 4, unknown keyword

### 9.6. service (Print Service Information to Terminal)

Syntax:            ?service or service  
Parameter:        none

Description:       The instruction service returns a multi-line parameter and state list of the currently used controller settings. It may be used for debugging or in case of service requests with a terminal program.

Response:         Many lines of text including e.g. serial number, parameters, states etc.

Example:           service

## 9.7. pci (Is PCI Bus)

Syntax:           ?pci or pci

Parameter:       none

Description:      The instruction pci returns:

0 = Controller is a desktop version

1 = Controller is a PCI card and plugged in a PCI slot

Response:        0 or 1

Example:          pci => 0

## 10. General Adjustments

With the following instructions the parameters of the controller are widely scalable to the given mechanic construction and to customer requirements. The controller is adaptable to the requested requirements.

### 10.1. dim (Unit for Positions)

Syntax:                   !dim or ?dim  
Parameter:               x, y, z or a  
                            0 to 8

Description:             The dim instruction sets the unit (or "dimension") of all input and output parameters related to length, e.g. position or move commands.

The provided units for length (parameters for dim) are:

0 => Micro steps  
1 =>  $\mu\text{m}$   
2 => mm (Tango default)  
3 =>  $360^\circ$   
4 => revolutions  
5 => cm  
6 => m  
7 => inch  
8 => mil

Examples:

```
!dim 4 1     the selected dimension for X is [revolutions] and for Y is [ $\mu\text{m}$ ].  
!dim z 2    the selected dimension for Z is [mm]  
?dim        responds the dimensions for all axes.  
?dim a      responds the dimension of the a-axis.  
dim 2 2 2 2 set dimension for all axes to [mm]
```

Response:     Current settings

Hint:         For dimensions 3 (=360°) and 4 (=revolutions) you should use a spindle pitch of 1mm to prevent rounding error.

### 10.2. pitch (Spindle Pitch)

Syntax:                   !pitch or ?pitch  
Parameter:                x, y, z or a  
                            0.0001 to 68

Description:             This instruction sends the spindle pitch (given by mechanic components) to the controller. It will be taken for all further calculations.

Response:                current spindle pitch

Examples:

```
!pitch 4.0 1.0    set spindle pitch X=4[mm] and Y=1[mm]  
!pitch z 2.0     set spindle pitch Z=2[mm]  
?pitch           query all axes for their spindle pitch  
?pitch a         query spindle pitch for a-axis
```

### 10.3. gear (Gear Ratio)

Syntax: !gear or ?gear  
Parameter: x, y, z or a  
0.001 to 1000

Description: This instruction transmits the gear ratio to and from the controller. The ratio is 1, if the motor is directly mounted on the spindle.

Response: current gear ratio

Examples:  
!gear 10 set gear ratio X=1/10  
!gear 4.0 1.0 set gear ratio X=1/4 and Y=1/1  
!gear z 10.0 set gear ratio Z=1/10  
?gear query all axes for their gear ratio  
?gear a query A-axis for its gear ratio

### 10.4. motorsteps (Motor Steps Per Revolution)

Syntax: !motorsteps or ?motorsteps  
Parameter: x, y, z or a  
[multiples of 4]

Description: This instruction sets or reads back the steps per revolution of the attached motor. Commonly the most motors have 200 steps per revolution (which is 1.8° full step). Other motors may have 400 or 500 steps per revolution. The motor steps parameter must be a multiple of 4.

Response: Currently used motorsteps

Examples:  
!motorsteps 200 200 400 set motor steps for X and Y to 200 and Z to 400  
!motorsteps x 500 set motor steps for X to 500  
?motorsteps query all axes for their motorsteps  
?motorsteps a query A-axis for its motorsteps

### 10.5. accel (Maximum Acceleration)

Syntax: !accel or ?accel  
Parameter: x, y, z or a  
0.01 to 20.00 [m/s<sup>2</sup>]

Description: This instruction transmits the maximum acceleration, which is currently used or shall be used for calculation of all consecutive moves.

Response: current acceleration in m/s<sup>2</sup>

Examples:  
!accel 1.00 set acceleration X=1.0[m/s<sup>2</sup>]. Other axis are not affected.  
!accel 1.00 1.50 set acceleration X=1.0[m/s<sup>2</sup>] and X=1.5[m/s<sup>2</sup>]  
!accel z 0.2 set acceleration Z=200[mm/s<sup>2</sup>]. Other axis are not affected.  
?accel query all axes for their current acceleration.  
?accel z query Z axis for its acceleration.

## 10.6. vel (Velocity)

Syntax: !vel or ?vel  
Parameter: x, y, z or a  
0.0000025 to 100

Description: This instruction transfers the velocity in motor revolutions per second, which is currently used or shall be used for calculation of all consecutive automatic moves.

Response: Currently selected velocity

Examples:

```
!vel 10          set velocity X=10[revolution/s]. Other axis are not affected.
!vel 1.0 15      set velocity X=1[revolution/s] and Y=15[revolution/s].
!vel z 0.1       set velocity Z=0.1[revolution/s].
?vel            query all axes for their velocities.
?vel x          query x axis for its velocity.
```

Additional information: The maximum velocity for automatic point to point moves has a resolution of 0.0001[revolution/s]. A query has a resolution of 0.001[revolution/s]. Nevertheless the internal representation and calculation is done with the higher resolution of the commanded number. If you need the higher accuracy also for queries, we recommend to use the command velfac in addition.

## 10.7. velfac (Velocity Factor)

Syntax: !velfac or ?velfac  
Parameter: x, y, z or a  
0.01 to 1.00

Description: This instruction transfers the velocity factor, which is currently used or shall be used for calculation of all consecutive automatic moves. It is internally multiplied to the velocity (vel).

Response: Currently used velocity factor [0.01 to 1.00]

Examples:

```
?velfac z        query X axis for its current velocity factor.
!velfac x 0.1    set velocity X to 1/10 of current velocity.
?velfac         query all axes for their current velocity factors.
```

## 10.8. secvel (Secure Velocity)

Syntax: !secvel or ?secvel  
Parameter: x, y, z or a  
1 to 100 [mm/s]

Description: The security speed limitation is used as long as the axis is not calibrated and range measured ("cal", "rm"). The velocity unit is always mm/s and does not depend on the "dim" state. It prevents the microscope stage from mechanical damage as long as the controller does not know the mechanical limits. It may also be used as a workaround, if running a cal/rm is not wanted.

Response: Currently used secure velocity [1 to 100 mm/s]

Examples: !secvel 100 100 100 => Set maximum possible velocity of X Y Z

## 10.9. maxcur (Query Maximum Motor Current)

Syntax: ?maxcur

Parameter: x, y, z or a

Description: This instruction reads the maximum possible motor current.

Response: maximum motor current in Ampere [A]

Examples:

?maxcur y query Y axis for its maximum motor current  
 ?maxcur query all axes for their maximum motor currents

## 10.10. cur (Motor Current)

Syntax: !cur or ?cur

Parameter: x, y, z or a  
 0.2 to [maximum current]

Description: This instruction transfers the current used or required motor current. The maximum current is limited by hardware and may be checked by the "maxcur" command.

Response: Selected motor current in Ampere

Examples:

!cur 1.0 set X motor current to 1[A]  
 !cur 1.0 2 set motor current for X=1[A] and Y=2[A]. Other axes are not affected.  
 !cur z 0.3 set Z motor current to 0.3[A].  
 ?cur query all axes for their motor currents.  
 ?cur x query X for its motor current.

## 10.11. reduction (Motor Current Reduction Factor)

Syntax: !reduction or ?reduction

Parameter: x, y, z or a  
 0 to 1.0

Description: This instruction transfers the current reduction factor. When the axis is idle, the motor current is reduced by this factor. A value of 1 disables the reduction, so even when the axis is stopped the motor current remains the same. Motor current reduction can be used to keep the motor temperature low, but as a side effect it may add noise, decrease performance and position accuracy. Please note that when setting the reduction back to 1, this will first take effect after the next move. Please also refer to the "curdelay" command.

Response: Selected reduction(s) [0.00 to 1.00]

Examples:

!reduction .1 .7 Set idle currents X=0.1\*cur[A] and Y=0.7\*cur[A]  
 !reduction z 0.5 Set Z idle current to 0.5\*cur[A]  
 ?reduction Query all axes for their current reduction factors.  
 ?reduction x Query X for its reduction factor.



## 10.12. curdelay (Delay for Current Reduction)

Syntax:                   !curdelay or ?curdelay  
Parameter:               x, y, z or a  
                          0 to 10000 [ms]

Description:             At the end of each move the axis enters the idle state. If the motor current reduction factor is set to a value less than 1.0 this reduction will take effect after the curdelay time.

Response:                Selected delay of current reduction in [ms]

Examples:

```
!curdelay 100 300 set delay for motor current reduction X=100[ms] and Y=300[ms]
!curdelay z 450   set delay for motor current reduction Z=450[ms]
?curdelay        query all axes for their motor current reduction delay.
?curdelay x      query X-axis for its motor current reduction delay.
```

## 10.13. axis (Enable, Disable, Switch Off Axis)

Syntax:                   !axis or ?axis  
Parameter:                x, y, z or a  
                          -1, 0, 1

Description:             This instruction enables, disables, switches off axes. Or reads its current state. A disabled axis still powers the motor with its current, while a switched off axis loses its torque.

Response:                Current axis state (1=enabled, 0=disabled, -1=power stage off)

Examples:

```
!axis 1 1 1 1     enable all axes.
?axis x           query X-axis for its state.
!axis 1 0 1 0     disable Y and A axis while X and Z are enabled.
!axis y -1        switch off Y axis: power stage Y off.
?axis            query all axes for their state.
```

## 10.14. axisdir (Axis Direction)

Syntax: !axisdir or ?axisdir  
 Parameter: x, y, z or a  
 0 or 1

Description: This command reverses the specified axes. The meaning of the limit switches (E0 and EE) is also automatically exchanged against each other:

The E0 and EE switches are treated as virtual switches, E0 is the switch in negative direction, EE in positive. The hardware is reassigned to the opposite switch. Also for: swact, swpol, swtyp, readsw. Exception: The 'swin' function is not affected. Please make sure to first set the desired axis direction before setting the end switch types, polarity etc.!

It is not recommended to change direction during operation!

0 = Normal direrction, CAL switch => E0, RM switch => EE  
 1 = Reversed direrction, CAL switch => EE, RM switch => E0

Response: Current axis direction is 0=not changed or 1=changed

Examples:

!axisdir 0 1 0 1 Axis directions of Y and A are reversed.  
 ?axisdir x Query X for its current axis direction.

## 10.15. stopaccel (Emergency Stop Deceleration)

Syntax: !stopaccel or ?stopaccel  
 Parameter: 0.01 to 20 m/s<sup>2</sup>

Description: This instruction sets the deceleration for emergency stop conditions. It will be used by abort commands or when detecting an unexpected limit switch (e.g. no cal/rm move was performed).

Response: Currently used deceleration for stop

Example: !stopaccel 0.8 set the X stop deceleration to 0.8[m/s<sup>2</sup>]

## 10.16. motortable (Motor Correction Table)

Syntax: !motortable or ?motortable  
 Parameter: x, y, z or a  
 0 or number specified by factory

Description: This instruction adds a motor correction. The motor has to be measured for the specific application by factory. There a table number will be assigned and the customer may activate it by setting the corresponding motortable number. Using a wrong motortable will lead to increased noise and position error.

0 = No correction

Response: Currently used motortable(s)

Example:

!motortable 1 1 2 0 Select motortable 1 for X and Y, 2 for Z and no for A  
 !motortable x 0 Disable correction in x  
 ?motortable Returns the currently used tables for all axes

## 11. Limit Switch Instructions (Hardware and Software)

### 11.1. lim (Software Limits)

Syntax:           !lim or ?lim  
Parameter:       x, y, z or a  
                  + maximum position range

Description:       This instruction sets the maximum allowed positioning range. The upper and lower software limits shall be send together in a single !lim instruction. Remember: The unit (dimension) of the transmitted numbers depends on the value of instruction dim.

Response:         Currently used software limits

Examples:  
!lim 1000 1000 2000 2000       set the software limits for X and Y.  
!lim z -500 1700               set the software limits for Z.  
?lim                             query all axes for their limits.  
?lim a                           query A-axis for its limits.

### 11.2. limctr (Enable or Disable Limit Control)

Syntax:           !limctr or ?limctr  
Parameter:       x, y, z or a  
                  0 or 1

Description:       This instruction enables or disables the limit control or shows the current state of limit control. Attention, be careful: If limit controls are disabled, the controller doesn't care about limits. In this case the controller may damage system components. Limit control is enabled by default from power on.

Response:         Limit control state (0 = not active, 1 = active)

Example:  
!limctr y 0                     disable Y limit control, Y axis limit switches are ignored  
!limctr 1 1 1                  enable X,Y and Z limit control  
!limctr z 1                     enable Z limit control  
?limctr a                       query A-axis for its status of limit control  
?limctr                         query all axes for their status of limit control

### 11.3. nosetlimit (Do not set limits by cal/rm)

Syntax:               !nosetlimit or ?nosetlimit  
Parameter:           x, y, z or a  
                      0 or 1

Description:         This command enables or disables the setting of software limit switches during calibration and range measure. The default is nosetlimit=0 which means that the software limits are set by the cal/rm moves to these min/max positions.

Response:  
0 = set software limits to !cal and !rm positions  
1 = do not change software limits after !cal and/or !rm

Examples:  
!nosetlimit 1 1     X and Y axis do not take software limits after !cal and !rm  
!nosetlimit y 1    Y axis is does not set software limits of !cal and !rm move  
?nosetlimit        query all axes for their nosetlimit state  
?nosetlimit a      query A axis for its nosetlimit state

### 11.4. swtyp (Type of Limit Switch)

Syntax:               !swtyp or ?swtyp  
Parameter:           x, y, z or a  
                      0 or 1

Description:         This instruction transmits the type of the limit switches. The sequence is E0 REF EE for all axes. The REF switch currently not used by the Tango controller. Important: When using no axis parameter (x,y,z or a), the 3 values will be used for all axes! To set individual axes, please do this separately, use the axis parameter x,y,z or a. Please note that the E0 and EE switch are reassigned by the 'axisdir' command.

0 = PNP, which adds a pull-down resistor to the switch input  
1 = NPN, which adds a pull-up resistor (default)

Response:            Currently selected type

Examples:  
!swtyp 1 0 1        Set all limit switches to NPN type  
!swtyp z 0 0 1      Set Z axis limit switches E0=PNP, REF=don't care, EE=NPN  
?swtyp y            Query Y axis for its switch type

## 11.5. swpol (Polarity of Limit Switch)

Syntax:               !swpol or ?swpol  
 Parameter:           x, y, z or a  
                       0 or 1

Description:        This instruction transmits the polarity of the limit switches. The sequence is E0 REF EE for all axes. The REF switch currently not used by the Tango controller. Important: When using no axis parameter (x,y,z or a), the 3 values will be used for all axes! To set individual axes, please do this separately, use the axis parameter x,y,z or a. Please note that the E0 and EE switch are reassigned by the 'axisdir' command.

0 = switch has active low signal  
 1 = switch has active high signal

Response:           current polarity of the limit switches

Examples:

```
!swpol y 1 1 1     set polarity of Y limit switches (E0 REF EE) to positive edge.
!swpol 1 0 1      set polarity of limit switches (E0 REF EE) for all axes.
!swpol z 0 0 0     set polarity of Z limit switches (E0 REF EE) to negative edge.
?swpol a           query limit switch polarity of the A axis
```

## 11.6. swact (enable or disable limit switches)

Syntax:               !swact or ?swact  
 Parameter:           x, y, z or a  
                       0 or 1

Description:        This command enables or disables the limit switches. The sequence is always:

E0 REF EE

0 = switch is inactive (actuation state is ignored)  
 1 = switch is active

The REF switch is not used by the Tango controller. Disabling limit switches may damage your system. When using no axis parameter, the 3 values will be used for all axes! To set individual axes please do this separately, use the axis parameter x, y, z or a. If a switch is set to inactive, it is not necessary to cal or rm this switch to abolish the secvel speed limitation. Inactive switches always return a non actuated state when using the '?readsw' command. But the 'swin' command still returns the switches TTL logic level state. Please note that the E0 and EE switch are reassigned by the 'axisdir' command.

Response:           current state, if limit switches are enabled or disabled

Examples:

```
!swact 1 0 1      Enable cal and rm limit switches for all axes (REF disabled)
!swact z 1 0 1    Eet Z limit switches E0=enabled REF=disabled EE=enabled
?swact a          Query A-axis, if limit switches are enabled or disabled.
```

## 11.7. readsw (Read Status of Limit Switches)

Syntax: ?readsw  
Parameter: none

Description: This instruction reads the limit switch state of all axes.  
The response is a string of 12 characters, either 0 or 1.

0 = limit switch is currently not actuated or disabled  
1 = limit switch is currently actuated (axis is in switch)

In opposite to the "swin" command, readsw returns the active/inactive state and not the signal level of the switch. Also the readsw exchanges E0 and EE switches when axisdir is changed.

Please note that the switch state is only valid when the swtyp, swpol parameters are set correctly and the switch is activated by swact.

Sequence of the 12 characters is:

Axis: x y z a x y z a x y z a  
Switch: E0 E0 E0 E0 Ref Ref Ref Ref EE EE EE EE

E0 = lower limit switch (!cal command)  
Ref = Reference switch  
EE = upper limit switch (!rm command)

Response: Actuation state of limit switches, 0 if disabled

Examples: ?readsw => query all limit switch actuation states

## 11.8. swin (Read Limit Switch Input Level)

Syntax: ?swin or swin  
Parameter: none

Description: This instruction reads the limit switch signal directly.  
The response is a string of 12 characters, either 0 or 1.

0 = limit switch input signal is TTL low  
1 = limit switch input signal is TTL high

In opposite to the "readsw" command, swin reflects the TTL input levels. Also disabled switches are represented with their current TTL input signal level. Swin is not affected by the axisdir command (does not exchange E0 and EE switches). The Ref signals are not used.

Sequence of the 12 characters is:

Axis: x y z a x y z a x y z a  
Switch: E0 E0 E0 E0 Ref Ref Ref Ref EE EE EE EE

E0 = lower limit switch (!cal)  
Ref = Reference switch (!ref)  
EE = upper limit switch (!rm)

Response: Switch input signal TTL levelstate of limit switches

Examples: swin => query all limit switch signal levels

## 11.9. statuslimit (Limit Status)

Syntax: ?statuslimit or statuslimit

Parameter: none

Description: The instruction statuslimit responds a string with 16 characters. They represent the state of the software limits.

They are arranged in 3 groups:

- 0 - 3: Group 1 => cal state of axis 0-3 (x,y,z,a)
- 4 - 7: Group 2 => rm state of axis 0-3 (x,y,z,a)
- 8 - 11: Group 3 => lower software limit state of axis 0-3 (x,y,z,a)
- 12 - 15: Group 4 => upper software limit state of axis 0-3 (x,y,z,a)

The characters may have one of four values:

- => the software limit has not yet been modified since power on
- A => axis is already calibrated (!cal)
- D => axis is already range measured (!rm)
- L => software limit has been modified by (!lim)

Example: Assume ?statuslimit returns the string "AAA-DLLLL"

This means in detail:

```
[ 0] A -> X-axis is calibrated
[ 1] A -> Y-axis is calibrated
[ 2] - -> Z-axis is not calibrated
[ 3] A -> A-axis is calibrated
[ 4] - -> X-axis is not range measured
[ 5] - -> Y-axis is not range measured
[ 6] - -> Z-axis is not range measured
[ 7] D -> A-axis is range measured
[ 8] - -> X-axis lower software limit is not modified
[ 9] L -> Y-axis lower software limit is modified
[10] L -> Z-axis lower software limit is modified
[11] - -> A-axis lower software limit is not modified
[12] L -> X-axis upper software limit is modified
[13] - -> Y-axis upper software limit is not modified
[14] - -> Z-axis upper software limit is not modified
[15] L -> A-axis upper software limit is modified
```

## 12. Calibration and Range Measure Instructions

After each power on or '!reset' of the controller the operator may run a calibration (instruction !cal) followed by a range measure (instruction !rm), if the system is equipped with the corresponding limit switches.

This also disables the travel speed limit set by 'sevel'. So the controller is able to smoothly stop the axes automatically before they crash into the microscope stage hardware limits.

These instructions set the limits very close to the limit switches. An additional offset for the these limits may be specified with the instructions !caliboffset and !rmoffset. Depending on the end switch version (e.g. light barrier), adding a position offset of 1 up to 10 millimeter may lead to a more stable !cal / !rm result.

### 12.1. cal (Command a Calibration)

Syntax: !cal or cal  
Parameter: x, y, z or a

Description: This instruction moves either the specified or all currently enabled axes in negative direction towards lower positions, until the limitswitch E0 is detected. It then moves with !calbspeed in plus direction out of the switch. If caloffset=0, the axis will stop moving as soon as the limitswitch E0 is not longer active, and set the reached position to 0. If caliboffset>0, the axis will continue moving until this distance is reached, and then set the position to 0. In both cases the reached position will also be taken as lower software limit, if nosetlimit is 1.

Response: each commanded (and enabled) axis answers either 'A' after a successful calibration or 'E' if an error occurred.

Examples:

```
!cal          do a calibration for all enabled axes
cal y        do a calibration for Y axis only (if enabled)
```

### 12.2. rm (Command a Range Measure)

Syntax: !rm or rm  
Parameter: x, y, z or a

Description: This instruction moves either the specified or all currently enabled axes in positive direction towards higher positions, until the limitswitch EE is detected. It then moves with !calbspeed in minus direction towards lower positions. If rmoffset=0, the axis will stop moving, as soon as the limit switch EE is not longer active. If rmoffset>0, the axis will continue moving until the distance defined by rmoffset is reached. In both cases the reached position will also be taken as lower software limit, if nosetlimit is 1.

Response: each commanded (and enabled) axis answers either 'D' after a successful range measure or 'E' if an error occurred.

Examples:

```
!rm          do a range measure for all enabled axes
rm z        do a range measure for Z axis only (if enabled)
```



### 12.3. caliboffset (Calibration Offset)

Syntax: !caliboffset or ?caliboffset  
 Parameter: x, y, z or a  
 Position

Description: This instruction specifies an extra offset position above the limitswitch E0 (towards higher positions) where to zero the axis and take this position as lower software limit. The unit depends on the current value of instruction dim. Allowed range is 0 to 30mm equivalent.

Response: Current calibration offset

Examples:

?caliboffset y query the Y-axis for its current calibration offset  
 !caliboffset 1 1 1 set the calibration offset to 1 for X,Y and Z.

### 12.4. roffset (Range Measure Position Offset)

Syntax: !roffset or ?roffset  
 Parameter: x, y, z or a  
 Position

Description: This instruction specifies an extra offset position below the limitswitch EE (towards lower positions) where to take the upper software limit. The unit depends on the current value of instruction dim. Allowed range is 0 to 30mm equivalent.

Response: current calibration offset

Examples:

?roffset z query the Z-axis for its current range measure offset  
 !roffset 1 1 1 set the range measure offset to 1 for X,Y and Z.

### 12.5. caldir (Calibration Direction)

Syntax: !caldir or ?caldir  
 Parameter: x, y, z or a  
 0 or 1

Description: This instruction set the calibration direction to either positive or negative positions. Default is negative direction. If set to positive(=1), the upper software limit is set. This instruction is not possible for systems with encoders.

Response: 0 = cal move to negative direction  
 1 = cal move to positive direction

Examples:

!caldir y 1 set Y axis calibration direction to positive  
 !caldir 0 0 1 set Z axis calibration direction to negative  
 ?caldir query all axes for their current calibration directions

## 12.6. calbspeed (Calibration Speed for Retraction)

Syntax: !calbspeed or ?calbspeed  
Parameter: range 1 to 100 [\*0.01 revolution/s]

Description: This instruction transmits the current or requested cal/rm calibration speed, which is taken for traveling out of the limit switch E0 and EE. There is only one value for all axes.

Response: current selected calibration speed

Examples:  
!calbspeed 15 set the retraction speed to 0.15 [revolutions/s] for all axes.  
?calbspeed query the controller for current retraction speed.

## 12.7. calrefspeed (Reference Signal Calibration Speed)

Syntax: !calrefspeed or ?calrefspeed  
Parameter: range 1 100 [\*0.01 revolution/s]

Description: This instruction transmits the current or requested reference calibration speed. This speed is taken for systems with encoders, when searching the reference on the scale. The default is 32. There is only one value for all axes. The value is not stored with !save instruction.

Response: current selected calrefspeed

Examples:  
!calrefspeed 5 set the retraction speed to 0.05 [revolutions/s] for all axes.

## 12.8. calpos (Calibration Position)

Syntax: !calpos or ?calpos  
Parameter: x, y, z or a  
position value

Description: This instruction is used for systems with encoders only. During calibration the encoder signal period is stored as soon as the E0 limit switch is left and may be queried later on with ?calpos.

This position may also be set to an other value. The value depends on the unit set by "dim". Allowed range is 0 to 30mm equivalent.

Response: within range of one encoder signal period

Examples:  
?calpos y query Y-axis for its calibration position  
!calpos 0 0 0 set calibration position to zero (X,Y and Z)  
?calpos query all axes for their read calibration position

## 12.9. refdir (Direction for Searching Reference Signal)

Syntax: !refdir or ?refdir

Parameter: x, y, z or a  
0 or 1

Description: This instruction is intended for systems with encoders and transmits the current or requested direction for searching the reference point on the scale. The default is 0 for minus direction.

Response: 0 = negative direction  
1 = positive direction

Examples:

!refdir y 1 set the Y-axis reference direction to positive direction  
?refdir query all axes for their reference directions

## 13. Move Instructions

All move instructions include an automatic linear interpolation. Axis, which are started together are reaching the destination at the same time. Nevertheless, user requested parameters, as specified for a single axis, are always calculated and the parameter given rules are not hurt. The user is free to also start the axis independent from each other. In this case each axis drives which its own maximum parameters and the axis do not reach the destination at the same time.

### 13.1. moa (Move Absolute)

Syntax:               !moa or moa  
Parameter:           x, y, z or a

Description:         This instruction moves one or more axes to a requested destination. The unit of the input numbers depends on instruction dim.

Response:            each commanded (and enabled) axis answers either '@' after an successful move or 'E' if an error occurred.

Examples:  
moa x 10.2            the X-axis is moved to position 10.2 ([mm] assume dim=2)  
moa 10 0 20          the axis X,Y,Z are moved to the given position  
moa y 34.5            the Y-axis is moved to position 34.5 ([mm] assume dim=2)

### 13.2. mor (Move Relative)

Syntax:               !mor or mor  
Parameter:           x, y, z or a

Description:         This instruction moves one or more axes relative to the current position. The unit of the input numbers depends on instruction dim.

Response:            each commanded (and enabled) axis answers either '@' after an successful move or 'E' if an error occurred.

Examples:  
mor x 12.3            the X-axis is moved by 12.3 ([ $\mu$ m] assume dim=1)  
!mor 1 1              the X and Y-axis are moved by 1 ([mm] assume dim=2)  
  
!mor a 298            the A-axis is moved by 298 (unit depends on dim)  
!mor -10 0 0 -10     the X and A-axis are moved by -10 (unit depends on dim)

### 13.3. m (Move Relative Shortcut)

Syntax: !m or m  
Parameter: none

Description: The instruction is a shortcut (abbreviation) of mor. It is useful to speed up communication especially for consecutive identical vectors. The vector is taken from the preceding !mor or !distance instruction. The instruction will move enabled axes if their distance is not zero.

Response: depends on state of autostatus.

This is one example of consecutive moves:

```
!moa 1 2 3 4    will position to 1 2 3 4
!mor 1 1 1 1    will move to 2 3 4 5
m              will move to 3 4 5 6
!distance 0 2 0 0
m              will move to 3 6 5 6
```

### 13.4. distance (Distance for m)

Syntax: !distance or ?distance  
Parameter: x, y, z or a  
Distance (+-2600mm max.)

Description: This instruction transmits the travel distance for !m instructions. The unit depends on the selected dimension (refer to "dim" instruction).  
Hint: The distance value is also set by executing a !mor command.

Response: current value for distance (unit depends on dim instruction)

Examples:

```
?distance      query current distance values for all axes
?distance z    query Z-axis for its distance value
!distance 10 20 set X and Y distance
!distance 1 2 3 set X,Y and Z distance
!distance y 20 set Y distance. Other axes keep their distance value.
```

### 13.5. moc (Move to Center)

Syntax: !moc or moc  
Parameter: x, y, z or a

Description: This instruction centers all enabled axis to the midpoint between lower and upper software limits. The recommended precondition is to first execute the instructions !cal and !rm.

Response: each successful centered axis responds with ,@'.

Examples:

```
moc          moves all axes to their centers
moc z       the Z-axis is centered, with no move of other axis
```

## 13.6. speed (Speed Move)

Syntax:           !speed or ?speed  
Parameter:       x, y, z or a  
                  + 100

Description:       This instruction moves one or more axes with requested speed, and not to a specified position. The command may be stopped by setting speed back to zero.  
                  As this is a "digital Joystick" function, this command only affects axes with Joystick enabled.

Response:         current value for speed [revolutions/s]

Examples:

```
!speed 33 0.01     set speed for X= 33[revolutions/s] and Y= 0.01[revolutions/s]
!speed 0           set speed for X-axis to 0[revolutions/s] (stop X axis)
!speed 10          set speed for X-axis to 10[revolutions/s]
!speed y 0.001     set speed for Y-axis
?speed             query all axes for their current speed
?speed z           query Z-axis for current speed
```

## 13.7. a (Abort the Current Move)

Syntax:           !a or a  
Parameter:        none

Description:       This instruction stops all axes and sets them into position reached state.  
                  You may also send a "Ctrl+C" (hex 0x03) command instead.

Response:         Each configured axis responds an ,@'.

Example:           a

## 13.8. delay (Set the Delay Time for Consecutive Moves)

Syntax:           ?delay or !delay  
Parameter:        0 to 10000 [ms]

Description:       This instruction transmits the current or requested delay time, which shall be kept before an automatic move starts.

Response:         Delay time currently used, in [ms]

Examples:

```
!delay 500         Set the delay time to 0.5 seconds
?delay             Query the current delay time
```

## 13.9. pause (Set the Pause after Position Reached)

Syntax: ?pause or !pause

Parameter: 0 to 10000 [ms]

Description: The pause time by which the "@@@" autostatus response is delayed. This may be used to insert an automatic settling time after a move command. It is the complement to the delay function.

Response: Pause time currently used, in [ms]

Examples:

!pause 10 Delay the autostatus response by 10 milliseconds

?pause Query the current pause time

## 13.10. pos (Read or Set Position)

Syntax: !pos or ?pos

Parameter: x, y, z or a  
Position (+- 2600mm max.)

Description: This instruction either reads or sets the current position. The used position unit depends on the selected dimension (refer to "dim" instruction). If an encoder and the 'encpos' is enabled, it returns the encoder position of the axis.

Response: Axis position(s) (depends on dim, and enc/encpos state)

Examples:

?pos Query all axes for their positions

!pos100 200 Set positions of X=100 and Y=200 (unit depends on dim)

!pos 0.1 Set the position X=0.1 (unit depends on dim)

!pos y 2000 Set the position Y=2000 (unit depends on dim)

?pos z Query Z-axis for its position

## 14. Joystick, Tackball and Handwheel Instructions

All manual moves with human input devices, like joystick and trackball, are limited by the limit switches automatically. You are allowed to unplug, plug and exchange these input devices, while the controller is switched on (hot plug). You can imagine the instruction speed as a digital, accurate joystick, nevertheless this instruction is an automatic controlled move. It is allowed to drive axis in different operation modes, e.g. to move X and Y axis with a moa instruction, while Z is controlled with speed instruction. This feature is basic and provides e.g. dynamic auto focus while moving the table.

### 14.1. joy (Generally Enable/Disable/Set Joystick Mode)

Syntax:                   !joy or ?joy  
 Parameter:               0, 1, 2, 3, 4 or 5

Description:  
 !joy 0                    disable the joystick/trackball mode  
 !joy 2                    enables joystick/trackball with position counting  
 Using other values than 0 or 2 is not recommended.

Response:                Current joystick mode

Examples:  
 !joy 2                    set joystick mode 2 (on)  
 ?joy                     query current mode

### 14.2. joydir (Joystick Direction or Assign Joystick)

Syntax:                   !joydir or ?joydir  
 Parameter:               x, y, z, or a  
 0, +1, +2,

Description:             This instruction sets the joystick/trackball direction or disables the joystick/trackball. Please note that when using a 4 axis controller with a 3 axis Joystick, its 3<sup>rd</sup> axis may be assigned to Z and/or A by enabling one or both Z, A joydirs. A value of 0 disables the joystick/trackball for this axis. A value of +-1 enables the joystick with full current. A value of +-2 enables the joystick with reduced current function support (refer to reduction/curdelay). Please make sure that the joystick function is globally enabled by the joy command.

Response:                current joydir settings

Examples:  
 !joydir 1                enable X-axis joystick reversed without current reduction.  
 ?joydir                 query all axes for their joystick direction settings.  
 !joydir 2 2 0 2         set positive direction, reduced current, assign the joysticks 3<sup>rd</sup> axis to the controller A axis only.



### 14.3. joywindow (Joystick Window)

Syntax: !joywindow or ?joywindow  
Parameter: 0 to 100

Description: This instruction sets the with of middle range, where joystick deflection has no effect to motor movement.  
Please note that there is only one value for all axes!  
This value should not be reduced, as this may result in slow unwanted moving of axes even when the joystick is apparently not deflected. Increasing the value will result in a loss of speed resolution.

Response: current value for joywindow

Examples:

?joywindow query for current joystick window  
!joywindow 14 set joystick window to 14

### 14.4. joyspeed (Joystick Speed Presets for BPZ Device)

Syntax: !joyspeed or ?joyspeed  
Parameter: 1, 2 or 3 and  
0.0001 to 100 [revolutions/s]

Description: This instruction sets the joystick speeds for the three BPZ device speed buttons only (Slow, Medium, Fast). Unit is in motor revolutions per second (as 'vel' instruction). Each button has to be set individually:

1 = Slow Button  
2 = Medium Button  
3 = Fast Button

Response: Current speed assigned to the button

Examples:

?joyspeed 1 Query for "Slow" joystick button speed  
!joyspeed 3 30 Set "fast" joystick button speed to 30 [revolutions/s]

### 14.5. key (Read HDI Device Key State)

Syntax: ?key or key  
Parameter: none or key number (1,2,3,4)

Description: This instruction reads the state of up to 4 HDI device keys.  
0 = key is currently released or not available  
1 = key is currently pressed

Response: 1 or 4 Key states, each either 0 or 1

Examples: key => query all keys, returns 4 numbers, e.g. 0 0 0 0  
key 1 => query only key 1 (e.g. F1 Joystick Key)

## 14.6. hwfactor (Handwheel Transmission Factor)

Syntax: !hwfactor or ?hwfactor  
Parameter: none or x, y, z, a  
and -200 to +200

Description: This instruction sets or reads the handwheel transmission factor, which is a floating point number between -200.0 and +200.0. A sign change may be used to change direction. The factor represents the stage travel distance in millimeter per handwheel knob revolution. Please note that when a higher transmission factor is selected, the step resolution becomes more coarse.

Response: Currently used handwheel factor(s)

Examples:

```
!hwfactor 10 10 => One knob revolution in X or Y results in 10mm travel
!hwfactor X 100 => One knob revolution in X results in 100mm travel
?hwfactor       => Query all axes for their transmission factor
```

## 14.7. hwfilter (Handwheel Noise Filter)

Syntax: !hwfilter or ?hwfilter  
Parameter: 0 or 1

Description: This instruction sets or reads the handwheel noise filter state.

1 = Noise filter is active (recommended, default)  
0 = Noise filter is deactivated (finer step resolution)

The filter can only be activated/deactivated for all axes. Disabling the filter may result in some inaccuracy between automatic moves, as the noise will cause slight position jitter when move commands are completed.

Response: Current state of handwheel filter

Examples:

```
!hwfilter 0 => No noise filter for handwheel, increased finer resolution
?hwfilter   => Query hwfilter state
```

## 14.8. tbfactor (Handwheel Transmission Factor)

Syntax: !tbfactor or ?tbfactor  
Parameter: none or x, y, z, a  
and -200 to +200

Description: This instruction sets or reads the trackball transmission factor, which is a floating point number between -200.0 and +200.0. A sign change may be used to change direction.

Response: Currently used trackball factor(s)

Examples:

```
!tbfactor X 100 => X axis is 10 times more sensitive than the default setting
?tbfactor       => Query all axes for their transmission factor
```

## 14.9. tvrjoy (Pulse and Direction Joystick Functionality)

Syntax: !tvrjoy or ?tvrjoy

Parameter: 0, z, a

Description: This instruction enables and assigns the AUX-IO pulse and direction input to an axis for simple joystick functionality. The behavior is similar to the trackball, which is available as HDI device.

Important: This option must not be used for absolute positioning of axes by an external controller. Please use the tvr functionality for this applications.

0 = Disabled  
z = Assigned to Z-axis  
a = Assigned to A-axis

Response: Currently assigned axis

Examples:

```
!tvrjoy 0          Disable AUX-IO tvr joystick function
!tvrjoy z          Assign AUX-IO tvr joystick function to Z-axis
?tvrjoy           Query assigned axis
```

## 14.10. tvrjoyf (Pulse and Direction Joystick Factor)

Syntax: !tvrjoyf or ?tvrjoyf

Parameter: -200 to +200

Description: This instruction sets or reads the tvrjoy transmission factor, which is a floating point number between -200.0 and +200.0. A sign change may be used to change direction.

Response: Currently used tvr factor

Examples:

```
!tvrjoyf 100       Axis is 10 times more sensitive than the default setting
?tvrjoyf           Query tvrjoy transmission factor
```

## 15. Digital and Analogue I/O

The Tango provides several digital I/O, two analogue outputs (channel 0 and 1) and one analogue input. These are available on the optional auxillary I/O port. The analogue output channel 2 is reserved for special purpose. Furthermore the HDI Interface analogue inputs may be read as well, if no HDI-device is connected.

### 15.1. digin (Digital Input)

Syntax:               ?digin or digin  
Parameter:           none or 0 to 15

Description:         This instruction queries the logic state of one or all digital inputs. If no parameter is used all inputs are returned as a string of 16 characters, ASCII 0 or 1, LSB (channel #0) first.

Response:            logic state of digital inputs

Examples:  
?digin                query all digital inputs  
?digin 8             query digital input 8

### 15.2. digout (Digital Output)

Syntax:               !digout oder ?digout  
Parameter:           string of 0s and 1s,  
                      0 to 15, 0 or 1

Description:         This instructions sets or reads back the logic level of the optional digital outputs. Outputs may be set either by a string of 0s and 1s or by channel number and signal level. The string is LSB first (channel 0 is the leftmost).

Response:            current output state

Examples:  
!digout 11110000    The digital outputs 0,1,2,3 are set to logic ,1' and the outputs 4,5,6,7 are set to logic ,0'.  
!digout 5 1         set digital output #5 to logic 1  
?digout             query the current state of all outputs  
?digout 8           query the current state of output 8  
!digout 7 0         set output 7 to 0

### 15.3. anain (Analogue Input)

Syntax: ?anain  
 Parameter: c (c = channel)  
 0 to 15 (channel number)

Description: This instruction reads the current value of one analogue input channel. The range is decimal from 0 (=0V) to 1023 (=5V).

Channel No	Connector	Pin	Signal Name
0	HDI	1	Joystick X
1	HDI	2	Joystick Y
2	HDI	3	Joystick Z
3	HDI	4	
4	HDI	5	Speedpoti
5	HDI	6	
6	HDI	7	
7	HDI	8	
8	HDI	9	
9	HDI	10	HDI-ID
10	AUX-IO	9	ANAIN0
11	internal	-	U-HIP
12	internal	-	V-MOT
13	EXT	20	X-ID0
14	EXT	18	X-ID1
15	internal	-	REF (2,5V)

Calculating the internal motor voltage:

$$U_{mot}[V] = (5 / 1023) * [anain\ c\ 12] * (55.7/4.7)$$

More accurate:

$$U_{mot}[V] = (2.5 / [anain\ c\ 15]) * [anain\ c\ 12] * (55.7/4.7)$$

Calculating the internal PSE voltage:

$$U_{mot}[V] = (5 / 1023) * [anain\ c\ 11] * (14.7/4.7)$$

More accurate:

$$U_{mot}[V] = (2.5 / [anain\ c\ 15]) * [anain\ c\ 11] * (14.7/4.7)$$

Calculating the case temperature (if available):

$$T[°C] = (250 / [anain\ c\ 11]) * [anain\ c\ 14]$$

Example:  
 ?anain c 10      Query level of channel 10 (analogue input of AUX-IO connector)

## 15.4. anaout (Analogue Output)

Syntax:           !anaout or ?anaout  
 Parameter:       0 to 100 in percent   (100% = 10V)  
                   c                           (c = single channel keyword)  
                   0, 1 or 2                 (single channel number)

Description:       This instruction sets and reads the values for analog outputs in percent. There are two ways to access the values, with or without the 'c' keyword (see examples below). So it is possible to address a single channel by using the 'c' or channel 0 or all channels by directly writing the percent values. Fractional numbers may be used, too. 100% corresponds to 10 Volts.

Channel No	Connector	Pin	Signal Name
0	AUX-IO	10	ANOUT0
1	AUX-IO	11	ANOUT1
2	reserved	-	-

Response:           Current value of analogue output in percent

Examples:  
 !anaout 100 50.1   Set channel 0 = 100% (10V) and channel 1 = 50.1% (5.01V)  
 !anaout 75         Set channel 0 = 75% (7.5V)  
 !anaout c 1 25.3   Set channel 1 to 25.3% (2.53V)  
 ?anaout            Query all channels for their output values  
 ?anaout c 0        Query channel 0 output value

## 15.5. stoppol (Polarity of Stop Input Signal)

Syntax:           !stoppol or ?stoppol  
 Parameter:       0 (= active low), 1 (= active high)

Response:           current polarity of stop signal

Description:       The stop input has an internal pull-up resistor to 5V. If you connect an NO (normal open) stop switch, you have to select low active. For NC (normal close) switches please select high active.

Example:  
 !stoppol 1         => Set the polarity of the AUX-I/O stop input to active high.

## 15.6. shutter (Shutter Out Signal of AUX-IO)

Syntax:           !shutter or ?shutter  
 Parameter:       0 (= low), 1 (= high)

Response:           Output level of shutter signal

Description:       Manually set the AUX-IO shutter out signal to the desired TTL level.

Example:  
 !shutter 1         => Set the shutter out signal into TTL high state.

## 16. Encoder Instructions

To enable encoder functionality, first the encoder mask has to be set for the corresponding axes. After that a cal move will activate the encoders ('enc'=1), so they can be used. Manually setting the encoders 'enc' state to 1 is not recommended. This may cause trouble with when in closed loop mode, and in case of analog encoders the signal correction will be missing also.

### 16.1. encmask (Encoder Mask)

Syntax: !encmask or ?encmask  
Parameter: x, y, z or a,  
0 or 1

Response: Encoder enable mask

Description: The instruction reads or sets the encoder globally enable mask. It is necessary to unmask encoders (=1) as a first step in order to use them later. This instruction does not activate the encoders, it just globally enables the usage of them. Please note: Encoders get detected and finally used after a successful calibration command 'cal'. Here the signal correction data is measured, too.

Example:  
!encmask 1 1 1 Globally enable encoders for X, Y and Z-axis  
!encmask z 0 Globally disable encoder for Z-axis  
?encmask Query encoder mask state for all axes

### 16.2. enc (Encoder Active)

Syntax: !enc or ?enc  
Parameter: x, y, z or a  
0 or 1

Response: Encoder active state

Description: This instruction may be used to determine if a 'cal' command has activated the encoders or not. In order to activate them, the encmask has to be set to 1 first. It is not recommended to activate encoders manually by setting enc to 1. Refer to the 'encmask' description for further information.

0 = Encoder is inactive (not used)  
1 = Encoder is activated

Example:  
?enc Globally enable encoders for X, Y and Z-axis  
?enc x Globally disable encoder for Z-axis  
!enc x 1 Manually activate the X-axis encoder (not recommended!)

### 16.3. encperiod (Encoder Signal Period)

Syntax: !encperiod or ?encperiod  
 Parameter: x, y, z or a  
 0.0001 to 1.000 [mm]

Response: Encoder signal period(s)

Description: This command reads or sets the encoder signal period. The unit is always [mm].

Example:

```
!encperiod 0.5 0.5 0.001    Set encoder period for X and Y to 500µm, Z to 1µm
!encperiod y 0.020         Set encoder period of Y-axis to 20µm
?encperiod                  Read encoder period of all axes
?encperiod x                Read encoder period of X-axis
```

### 16.4. encdir (Encoder Counting Direction)

Syntax: !encdir or !encdir  
 Parameter: x, y, z or a  
 0 or 1

Response: Encoder counting direction

Description: The encoder counting direction is set automatically by the calibration 'cal' move. These directions must not be changed afterwards.  
 Only if the axis should be used for e.g. relative measurement and not for closed loop, it is ok to change the direction manually.

0 = Encoder counting direction default  
 1 = Encoder counting direction reversed

Example:

```
!encdir 1 1 1              Reverse encoder counting direction for all axes
!encdir x 1                Reverse encoder counting direction for X-axis only
?encdir                    Query encoder counting direction for all axes
?encdir y                  Query encoder counting direction for Y-axis only
```

### 16.5. encctl (Encoder has TTL Signal)

Syntax: !encctl or ?encctl  
 Parameter: x, y, z or a  
 0 or 1

Response: Currently selected encoder signal type(s)

Description: This command reads or writes the currently selected type of encoder signal processing.

0 = Encoder has analog sin/cos signals  
 1 = Encoder has digital quadrature A/B signals (e.g. RS422)

Example:

```
!encctl 0 0 1             Y and Y axis encoders are analog, Z is digital A/B encoder
!encctl z 1               Set Z encoder signal processing to digital
?encctl                   Query all axes for their currently used signal type
?encctl x                 Query X-axis for its currently used signal type
```



## 16.6. encref (Use Encoder Reference Signal)

Syntax:               !encref or ?encref  
Parameter:            x, y, z or a  
                          0 or 1

Response:             Encoder reference signal utilized state

Description:          This functionality is currently not supported.  
  
                          0 = Encoder reference signal not utilized  
                          1 = Encoder reference signal utilized during calibration

Example:  
!encref 1 1 0         Utilize encoder reference signal for X and Y-axis  
!encref x 1           Utilize encoder reference signal for Y-axis  
?encref               Query Encoder reference signal utilized state of all axes  
?encref x             Query Encoder reference signal utilized state of X-axis

## 16.7. encnas (Use Encoder NAS Error Signal)

Syntax:               !encnas or ?encnas  
Parameter:            x, y, z or a  
                          0 or 1

Response:             Encoder NAS error signal usage state

Description:          Before enabling this functionality please make sure, that the  
                          connected encoder provides a NAS error signal.  
                          If enabled, a encoder NAS error also generates an internal  
                          'err' error state. The NAS input signals an encoder error  
                          state by a TTL low level.  
  
                          0 = NAS encoder input state is ignored (default)  
                          1 = NAS encoder input signal is used for extended error  
                              detection

Example:  
!encnas 1 1 0         Utilize encoder NAS signal for X and Y-axis  
!encnas x 1           Utilize encoder NAS signal for Y-axis  
?encnas               Query Encoder NAS signal utilized state of all axes  
?encnas x             Query Encoder NAS signal utilized state of X-axis

## 16.8. encnasstatus (Encoder NAS Error Signal State)

Syntax: ?encnasstate or encnasstate

Parameter: x, y, z or a

Response: Encoder NAS error signal state

Description: Before enabling this functionality please make sure, that the connected encoder provides a NAS error signal. If enabled, a encoder NAS error also generates an internal 'err' error state. The NAS input signals an encoder error state by a TTL low level.

0 = NAS capable encoder signals: no error

1 = NAS capable encoder signals: error

Example:

encnasstate Query NAS signal error state for all axes

encnasstate x Query NAS signal error state for X-axis only

## 16.9. encerr (Encoder Error State)

Syntax: !encerr or ?encerr

Parameter: x, y, z or a

0

Response: Encoder error state

Description: This command reads or resets the encoder error state. On error the encoder signal is invalid and a potentially running closed loop for the corresponding axis is switched off.

0 = No error, normal function

1 = Encoder error

Example:

!encerr 0 Reset encoder error

?encerr Read encoder error states of all axes

?encerr x Read encoder error states of X-axis only

## 16.10. encpos (Encoder Position)

Syntax: !encpos or ?encpos  
 Parameter: x, y, z or a  
 0 or 1

Response: Position output type

Description: If set to 1 and the encoder is activated (corresponding enc = 1), a '?pos' returns the encoder position. Refer to the 'pos' and 'enc' commands for further information.

0 = pos command reads the user position (default)  
 1 = pos command reads the encoder position (if encoder active)

Example:

```
!encpos 1 1 0    a 'pos' command returns the encoder position for X and Y-axis
                  (if encoders are active)
!encpos x 1      a 'pos' command returns the encoder position for the X -axis
                  (if encoder is active)
?encpos          Read position output type for all axes
?encpos x        Read position output type for X-axis only
```

## 16.11. hwcount (Hardware Counter)

Syntax: ?hwcount or hwcount  
 Parameter: x, y, z or a

Response: Hardware counter reading(s)

Description: Hwcount returns the position(s) of the independent TTL encoder counter. It is a digital counter that counts the signal slopes (4 per period) and does not provide signal interpolation. So one signal period corresponds to a counter reading of 4. See also the 'clearhwcount' command.

Example:

```
hwcount          Returns the counter readings of all axes
hwcount x        Returns the counter readings of X-axis only
```

## 16.12. clearhwcount (Clear Hardware Counter)

Syntax: !clearhwcount or clearhwcount  
 Parameter: x, y, z or a

Response: Reset hardware counter reading(s)

Description: This command resets the hardware counter(s) to zero.

Example:

```
clearhwcount     Reset hwcount position of all axes to zero
clearhwcount x   Reset hwcount position of X-axis to zero
```

## 17. MR Encoder Instructions

### 17.1. mra (MR Amplitude Correction Factor)

Syntax: !mra or ?mra  
 Parameter: x, y, z or a  
 0.8 to 1.2

Response: Currently used correction factor(s)

Description: This command reads or sets the cosine amplification correction factor of the analogue encoder signal (here: sin/cos amplitude ratio). This factor is calculated automatically on each calibration move 'cal' and should not be changed. If the axis is manually controlled and only used for relative measurement, so that no 'cal' is possible, the user may determine the ratio itself and then write it into mra for more accurate results. Please also refer to the mro command.

Example:  
 ?mra Read MR signal correction factor of all axes  
 ?mra x Read MR signal correction factor of X-axis only  
 !mra x 1.0095 Amplify the X cosine signal by \*1.0095 compared to the sine

### 17.2. mro (MR Amplitude Offset Correction Value)

Syntax: !mro or ?mro  
 Parameter: x, y, z or a  
 -2048 to +2048

Response: Currently used correction value(s)

Description: This command reads or sets the sine and/or cosine offset compensation value as 16bit signed digits. This factor is calculated automatically on each calibration move 'cal' and should not be changed. If the axis is manually controlled and only used for relative measurement, so that no 'cal' is possible, the user may determine the offset itself and then write it into mro for more accurate results. Please also refer to the mra command.

Example:  
 ?mro Read MR signal offset value sine and cosine for all axes  
 ?mro x Read MR signal offset value sine and cosine for X-axis only  
 !mro 48 -100 0 0 0 0 Set X offset to sin=48digit, cos=-100digit, Y, Z = 0  
 !mro y 16 -28 Set Y offset to sin=16digit, cos=-28digit  
 !mro y 16 Set only sine offset of Y encoder

### 17.3. mrp (MR Signal Peak-To-Peak Measuring Result)

Syntax: !mrp or ?mrp  
Parameter: x, y, z or a  
-2048 to +2048

Response: [sine max] [sine min] [cosine max] [cosine min] reult(s)

Description: This command reads or sets the sine and/or cosine peak values, measured since they were reset the last time.  
It is just a measurement and has no effect to the signal processing itself. The returned values are signed 16bit digits.

Example:

```
?mrp x Returns [x_sin max] [x_sin min] [x_cos max] [x_cos min]
?mrp Returns the above, but for all axes (up to 16 values)
!mrp x 0 0 0 0 Reset the peak-to-peak measurement for x
!mrp x 0 0 Reset only the X sine min, max values
!mrp 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 Reset measurement for all 4 axes
```

### 17.4. mrt (MR Signal Level)

Syntax: ?mrt  
Parameter: x, y, z or a  
1 to 32767

Response: [sine] [cosine] reult(s)

Description: This command reads the corrected sine and cosine A/D converter results as signed 16bit digits.  
If there is no count parameter transmitted, the command returns 10 measurement result lines per default.

Example:

```
?mrt x 1 Returns one line with [x_sin] [x_cos] signal digits
?mrt 1 Returns the above, but for all axes (up to 8 values):
[x_s] [x_c] [y_s] [y_c] [z_s] [z_c] if 3 axes are configured

?mrt y 2 Returns two lines with [y_sin] [y_cos] signal digits
?mrt Returns 10 lines with all axes (up to 8 values per line)
?mrt x Returns one line with [x_sin] [x_cos] signal digits
```

## 18. Closed Loop Instructions

The closed loop control positions the stage to the scale position. So the inaccuracy of the drive is compensated. The closed loop control circuit is activated by the "ctr" command. But first, in order to activate the encoders they have to be unmasked "encmask" and a calibration move "cal" has to be executed (which activates "enc" and detects the counting direction).

### 18.1. ctr (Control Enable)

Syntax:               !ctr or ?ctr  
Parameter:           x, y, z or a  
                      0,1,2,3,4

Description:         This instruction activates the closed loop circuit.  
                      0 = Closed Loop OFF  
                      1 = Closed Loop Auto OFF each time position is reached  
                      2 = Closed Loop always ON (currently not supported!)  
                      3 = (currently not supported!)  
                      4 = (currently not supported!)

Response:            Closed loop state(s)

Examples:  
!ctr 0 0 0 0         Closed loop off for all axes  
!ctr x 1            Closed loop for X-axis switches off after position reached  
?ctr                Query closed loop states of all axes  
?ctr x              Query closed loop state of X axis

### 18.2. ctrc (Control Call)

Syntax:               !ctrc or ?ctrc  
Parameter:            1 to 100 [ms]

Description:         This instruction reads or sets the controller call interval.  
                      Unit is milliseconds.

Response:            Closed loop control call interval in milliseconds.

Examples:  
!ctrc 1             Closed loop control is executed every millisecond  
?ctrc               Query closed loop call interval

### 18.3. ctrt (Control Timeout)

Syntax:               !ctrt or ?ctrt  
Parameter:            0 to 10000 [ms]

Description:         This instruction reads or sets the control timeout.  
                      This is the time the closed loop tries get to the desired  
                      scale position. Unit is milliseconds.

Response:            Closed loop control timeout in milliseconds.

Examples:  
!ctrt 3000         Closed loop controller tries to reach the position for 3ms  
?ctrt              Query closed loop timeout

## 18.4. twi (Target Window)

Syntax:               !twi or ?twi  
Parameter:            x, y, z, or a  
                          [value corresponding 0.0001 to 1 mm] in dim units

Description:         This instruction reads or sets the closed loop control target window width. While increasing this value leads to position variance, setting a too narrow window may result in oscillation and closed loop timeouts.  
                          The unit depends on 'dim'.

Response:            Closed loop target window.

### Examples:

!twi 0.001 0.001    Closed loop target window is 1 $\mu$ m (if dim=2) for X and Y-axis  
!twi y 0.005        Closed loop target window is 5 $\mu$ m (if dim=2) for Y-axis  
?twi                 Query all axes for their target window  
?twi z               Query Z-axis for its target window

## 19. Trigger Signal Configuration

These commands read or modify the parameters for the trigger output signal. It may be used for synchronization of an external device like e.g. a video camera. The trigger output signal is available on the optional AUX-I/O connector. Access permission to the trigger functionality has to be enabled by factory. Before enabling the trigger function (by “!trig 1”), please make sure that all trigger settings have been made.

```
Example1:  !trig 0[CR]           Disable trigger
           !trigm 0[CR]        Choose trigger mode 0
           !triga x[CR]        Choose X axis as trigger source
           !trigd 0.100[CR]    Set trigger distance to 100µm (if dim = 2)
           !trigs 400[CR]     Set trigger pulse width to 0.4ms
           !trig 1[CR]        Enable trigger, set start position
```

```
Example2:  !trig 0[CR]           Disable trigger
           !trigs 120[CR]       Set trigger pulse width to 120µs
           !trigf 2500[CR]     Set pulse frequency to 2.5kHz
           !trigm 100[CR]     Choose trigger mode 100 (periodic signal)
           !trig 1[CR]        Enable trigger, set start position
```

Optional the “trigcount 0” command may be executed to reset the event counter.

### 19.1. trig (Trigger)

```
Syntax:      !trig or ?trig
Parameter:   0 (= disabled) or 1 (= enabled)
```

Description: This instruction enables or disables the trigger circuit. “!trig 1” also sets the trigger start position.

```
0 = Trigger function globally disabled
1 = Trigger function globally enabled
```

```
Response:   0 or 1
```

```
Examples:
!trig 0           Disable trigger circuit
?trig            Query for current state of trigger circuit
```

### 19.2. triga (Trigger Axis)

```
Syntax:      !triga or ?triga
Parameter:   x, y, z or a
```

Description: This instruction selects the axis on which to trigger

```
Response:   x, y, z or a
```

```
Examples:
!triga x         Select X-axis as trigger source
!triga y         Select Y-axis as trigger source
?triga          Query current trigger axis
```



### 19.3. trigm (Trigger Mode)

Syntax: !trigm or ?trigm  
 Parameter: 0 to 11, 100 to 103

Description: This instruction selects the required trigger mode.

Trigger Mode	Trigger Generation	Trigger Signal	Remarks
0		High active	First pulse when move starts
1		High active	First pulse when move starts
2		High active	First pulse when move starts
3	-- See Mode 0 --	Low active	Same as 0, signal inverted
4	-- See Mode 1 --	Low active	Same as 1, signal inverted
5	-- See Mode 2 --	Low active	Same as 2, signal inverted
6		High active	Triggers shifted by trigd/2
7		High active	Triggers shifted by trigd/2
8		High active	Triggers shifted by trigd/2
9	-- See Mode 6 --	Low active	Same as 6, signal inverted
10	-- See Mode 7 --	Low active	Same as 7, signal inverted
11	-- See Mode 8 --	Low active	Same as 8, signal inverted
100	Generates periodic trigger signals with the frequency chosen by the "trigf" parameter.	High active	Does not depend on position
101	-- See Mode 100 --	Low active	Same as 100, signal inverted
102	Allows manual forced trigger signals by the "trigger" command.	High active	Does not depend on position or time
103	-- See Mode 102 --	Low active	Same as 102, signal inverted

Response: Trigger mode as integer: 0 to 11, 100 to 103  
Examples: !trigm 3 Set Trigger Mode 3  
?trigm Query current trigger mode

## 19.4. trigger (Force Trigger Signal)

Syntax: !trigger or trigger  
Parameter: None

Description: This instruction generates a trigger output pulse. It is available in trigger modes 102 and 103. The pulse width is depending on "trigs" value.

Response: None

Examples:  
trigger Force trigger pulse now  
!trigger The same as above

## 19.5. trigs (Trigger Signal Length)

Syntax: !trigs or ?trigs  
Parameter: 0 to 2621400 [ $\mu$ s] (as multiples of 40 [ $\mu$ s])

Description: This instruction is used to adjust the trigger pulse width from 40 microseconds to 2.6214 seconds in increments of 40. (0 = shortest trigger signal width)  
If the parameter is not a multiple of 40 it will be rounded down to the next lower multiple (e.g. 100 --> 80). When read back, the corrected value is returned (here: 80).

Response: 0 to 2621400 ( $\mu$ s), as multiple of 40

Examples:  
!trigs 40 Set Trigger pulse width to 40  $\mu$ s  
!trigs 2500000 Set Trigger pulse width to 2.5 s  
?trigs Query current trigger pulse width

## 19.6. trigd (Trigger Distance)

Syntax: !trigd or ?trigd  
Parameter: >0 to 5000000 (unit depends on dim of the selected axis)

Description: This instruction selects the required trigger distance. After passing an interval of trigd with the selected axis, a trigger signal is generated.

Response: Trigger distance

Examples:  
!trigd 3 Set trigger distance to 3mm (if dim of selected axis is 2)  
!trigd 0.010 Set trigger distance to 10 $\mu$ m (if dim of selected axis is 2)  
?trigd Query current trigger distance

## 19.7. trigf (Trigger Frequency)

Syntax:           !trigf or ?trigf  
Parameter:       0.01 to 12500

Description:      This instruction generates periodic trigger pulses at the desired frequency. It is available at trigger mode 100. Frequency resolution is 1/40 $\mu$ s.

Response:        Trigger frequency

Examples:  
!trigf 2500        Generate periodic trigger pulses with 2.5kHz (each 0.4ms)  
?trigf            Query current trigger frequency

## 19.8. trigcount (Trigger Counter)

Syntax:           !trigcount or ?trigcount  
Parameter:        0 to 2147483647

Description:      This instruction reads or manipulates the counted trigger events.

Response:        Number of executed triggers

Examples:  
?trigcount        Query current trigger count  
!trigcount 0      Clear trigger counter

## 20. Snapshot Signal Configuration

These commands read or modify the parameters for the snapshot input signal, which may be generated from an external device for synchronization purpose. The snapshot input signal is in the moment only available on Joystick button "F2". If snapshot is enabled, each snapshot event will store the current axis positions in an array. Access permission to the snapshot functionality has to be enabled by factory. Please globally enable the snapshot function first ("sns 1") after all snapshot settings have been made.

**Example: Three snapshot positions are captured**

sns = 3  
 sns =

Element Index	Position X	Position Y	Position Z	Position A
1	1.0000	1.2345	1.2345	0
2	2.0000	1.2345	1.2345	0
3	3.0000	1.2345	1.2345	0
4	<i>invalid</i>	<i>invalid</i>	<i>invalid</i>	<i>invalid</i>
5	<i>invalid</i>	<i>invalid</i>	<i>invalid</i>	<i>invalid</i>

...                    ...                    ...                    ...                    ...

200	<i>invalid</i>	<i>invalid</i>	<i>invalid</i>	<i>invalid</i>
-----	----------------	----------------	----------------	----------------

In **sns = 0** the next snapshot will add a new row at index 4.  
 Same does the **!sns** command.  
**?sns** will return positions of row 3.  
**?sna x 2** will return 2.0000

### 20.1. sns (Snapshot)

Syntax:                    !sns or ?sns  
 Parameter:                0 (=disable) or 1 (=enable)

Description:              This instruction enables or disables the snapshot input.

Response:                Snapshot state

Examples:  
 !sns 0                    Disable snapshot  
 !sns 1                    Enable snapshot  
 ?sns                      Query state of snapshot circuit

### 20.2. snsl (Snapshot Level / Polarity)

Syntax:                    !snsl or ?snsl  
 Parameter:                0 (=active low) or 1 (active high)

Description:              This instruction sets the required snapshot signal polarity.

Response:                Currently used snapshot polarity

Examples:  
 !snsl 0                    Set snapshot input to active low  
 !snsl 1                    Set snapshot input to active high  
 ?snsl                      Query current snapshot input polarity

### 20.3. snsf (Snapshot Filter)

Syntax:           !snsf or ?snsf  
Parameter:       0 to 100 [ms]

Description:      This instruction reads or modifies the snapshot filter time, which is used to suppress glitches or spikes on noisy signals.

Response:         Currently used snapshot filter time

Examples:  
!snsf 0            Disable input filter  
!snsf 10           Set snapshot filter time to 10 ms  
?snsf             Query snapshot filter time

### 20.4. snsrm (Snapshot Mode)

Syntax:           !snsrm or ?snsrm  
Parameter:        0 (=capture), 1 (=move), 2 (=extended move)

Description:      This instruction reads or sets the snapshot mode (default=0).

0 = Capture positions with Joystick key "F2"  
1 = Move to Positions with Joystick key "F2"  
2 = Extended move:  
    F1: Step/move through position list forward (pointer+1)  
        (wraps around at the last element)  
    F2: Step/move through position list backward (pointer-1)  
        (wraps around at the first element)  
    F3: Move to start of list (first element)  
    F4: Moves to "prehome" position with "vel",  
        then to "home" position with "sevel"

Position capture and moves are always executed on all active axes.

Response:         Currently selected snapshot mode

Examples:  
!snsrm 0           Set snapshot mode to capture  
!snsrm 1           Set snapshot mode to move  
!snsrm 2           Set snapshot mode to extended move  
?snsrm             Query current snapshot mode

### 20.5. snsc (Snapshot Counter)

Syntax:           !snsc or ?snsc  
Parameter:        --

Description:      This instruction reads the snapshot counter, which shows the counted snapshots (= current snapshot array entries). This instruction may also be used to reset the counter to zero.

Response:         Current snapshot array entries (= number of snapshot events)

Example:  
?snsc             Query the number of detected snapshots.  
!snsc             Clear snapshot counter

## 20.6. snsp (Snapshot Position)

Syntax: !snsp or ?snsp  
 Parameter: x, y, z or a

Description: This instruction reads or writes the snapshot position. Writing positions appends them to the current position array. Reading positions returns the last captured position (last array element).

Remark: The position data unit depends on selected dimension 'dim'.

Response: Snapshot position value(s)

Examples:

```
!snsp 100 200    Append snapshot position for X and Y
!snsp 10 20 30   Append snapshot position for X, Y and Z axis
!snsp y 2000     Append snapshot position to Y = 2000
?snsp           Query all axes for their last captured snapshot positions
?snsp z        Query Z axis for its last captured snapshot position
```

## 20.7. snsas (Snapshot Array)

Syntax: !snsas or ?snsas  
 Parameter: x, y, z or a  
 and entry index from 1 to 200

Description: This instruction reads or writes to the snapshot position array, which may contain up to 200 elements. For reading, a valid element index may have a value of 1 to maximum the current snapshot counter value 'snsc'. For writing an index of snsc+1 may be used to append a position element to the array (snsc then gets updated by +1).

Remark: The position data unit depends on selected dimension 'dim'.

Response: Snapshot array position(s)

Examples:

```
?snsas 1        Query 1st snapshot entry for all axes positions
?snsas 33       Query 33rd snapshot entry for all axes positions
?snsas z 99     Query 99th snapshot entry for Z-axis position
?snsas x 199    Query 199th snapshot entry for X-axis position
!snsas 0        Clear the entire snapshot array
!snsas x 1 20.5 Set X position of first element to 20.5 (e.g. mm if dim=2)
!snsas 2 10 10 10 Set all axis positions of second array entry to 10
```

## 20.8. prehome (Snapshot PreHome Position)

Syntax: !prehome or ?prehome

Parameter: x, y, z or a

Description: This instruction sets the prehome position used by the snapshot extended move. The unit of the input position depends on instruction dim.  
See "snsm" 2 for more details.

Response: Position value(s)

Examples:

!prehome x 10.2 Set prehome position X-value to 10.2 (e.g. [mm] when dim=2)

!prehome 10 0 20 Set prehome position X,Y,Z

?prehome x Read currently used prehome X-position

?prehome Read currently used prehome positions of all axes

## 20.9. home (Snapshot Home Position)

Syntax: !home or ?home

Parameter: x, y, z or a

Description: This instruction sets the home position used by the snapshot extended move. The unit of the input position depends on instruction dim.  
See "snsm" 2 for more details.

Response: Position value(s)

Examples:

!home x 10.2 Set home position X-value to 10.2 (e.g. [mm] when dim=2)

!home 10 0 20 Set home position X,Y,Z

?home x Read currently used home X-position

?home Read currently used home positions of all axes

## 21. Scan Mode

### 21.1. scanmode (Switch to Scan Mode)

Syntax: !scanmode or ?scanmode

Parameter: 0 or 1

Description: This instruction switches the Tango controller into scan mode. In this mode the vector velocity is constant and can be set by 'scanvel'.

0 = normal operation

1 = scan mode

Response: Controller mode.

Examples:

```
!vel 10          set velocity X=10[revolution/s]. Other axis are not affected.
!vel 1.0 15     set velocity X=1[revolution/s] and Y=15[revolution/s].
!vel z 0.1      set velocity Z=0.1[revolution/s].
?vel           query all axes for their velocities.
?vel x         query x axis for its velocity.
```

### 21.2. scanvel (Scanmode Vector Velocity)

Syntax: !scanvel or ?scanvel

Parameter: 0.000001 to 1000 [mm/s]

Description: This instruction transfers the scanmode vector velocity in millimeter per second. As this is a vector mode there is only one velocity parameter.

Response: Currently selected velocity

Examples:

```
!scanvel 10     Set scanmode vector velocity to 10mm/s
?scanvel       Query scan mode velocity.
```



## 22. Document Revision History

No.	Revision	Date	Changes	Remarks
01	A	03. July 2007	New layout, improved and corrected descriptions, added new instructions, re-sorted instructions	Based on Tango firmware revision 1.26
02	B	09. July 2007	Added new instructions	
03	prelim. C	27. July 2007	twi example corrected	
04	C	03. Sept 2007	Added snapshot functions	